# NanoPhotometer **NP80/N60/N50/C40**

## User Manual

Version 2.0.2

Software Version 2.0.12465

## Declaration of conformity for the NanoPhotometer (C40/N50/N60/NP80)

This is to certify that the Implen NanoPhotometer conforms to the requirements of the following directives:

2014/35/EU	Low Voltage Equipment Safety Directive
2004/108/EC	EMC Directive
IEC 60529	Protection class IP20
2002/95/EC	Restrictions on the use of certain Hazardous Substances in Electrical and Electronic Equipment (ROHS)
2012/19/EU	EC Directive on Waste Electrical and Electronic Equipment (WEEE) 2003/108/EC & 2008/34/EC. By ensuring this product is disposed of correctly, you will help prevent potential negative consequences for the environment and human health, which could otherwise be caused by inappropriate waste handling of this product.
FCC 47 CFR Part15 §15.107 and §15.109	
EN 301 489-1 V1.9.2	Radio and ancillary equipment for portable use (portable equipment); EUT Operating frequency range: 2.4 – 2.4835 GHz
EN 300 328 V1.8.1	Electromagnetic compatibility and Radio spectrum Matters (ERM); Wideband transmission systems; Data transmission equipment operating in the 2,4 GHz ISM band and using wide band modulation techniques; Harmonized EN covering the essential requirements of article 3.2 of the R&TTE Directive
EN 301 489-17 V2.2.1	Electromagnetic compatibility and Radio Spectrum Matters (ERM)
IEC 62133 and UN38.3	Battery certification and transport test
Standards to which conformity is declared, where relevant, are as follows:	
IEC/EN 61010-1:2012	Safety requirements for electrical equipment for measurement, control, and laboratory use. General requirements.
EN61326-1:2013	Electromagnetic compatibility- generic emission standard electrical equipment for measurement, control, and laboratory use.

Signed:

Dr. Thomas Sahiri  
Managing Director  
Implen GmbH

## Contents

<b>1. NANOPHOTOMETER AT A GLANCE.....</b>	<b>7</b>
NANOPHOTOMETER OVERVIEW .....	7
NANOPHOTOMETER NP80 .....	8
NANOPHOTOMETER N60 .....	9
NANOPHOTOMETER N50 .....	10
NANOPHOTOMETER C40 .....	11
INSTRUMENT REAR PANEL (NP80/N60/N50/C40).....	12
INSTRUMENT BOTTOM VIEW (NP80/N60/N50/C40).....	12
ACCESSORIES .....	13
STANDARD ACCESSORIES.....	13
OPTIONAL ACCESSORIES.....	13
CONNECTIVITY.....	17
NANOPHOTOMETER SPECIFICATIONS.....	18
<b>2. GETTING STARTED.....</b>	<b>19</b>
SPECTROPHOTOMETER INSTALLATION .....	19
SAFETY INFORMATION .....	19
UNPACKING AND POSITIONING .....	20
SOFTWARE INSTALLATION.....	21
NPOS OVERVIEW.....	21
REQUIREMENTS AND COMPATIBILITY.....	21
INSTALLING SOFTWARE ON COMPUTER.....	22
INSTALLING NANOPHOTOMETER APP ON TABLET OR SMARTPHONE.....	23
FIRST STEPS AND CONFIGURATION WIZARD.....	24
PRINTER INSTALLATION .....	24
<b>3. NANOPHOTOMETER BASICS.....</b>	<b>25</b>
APPLICATIONS OVERVIEW .....	25
ICONS .....	27
BUTTONS.....	28
SIDE TAB BAR .....	29
DATA PROCESSING DIALOGS .....	31
BASIC OPERATION.....	34
NANOVOLUME MEASUREMENT BASICS (N50/N60/NP80).....	34
CUVETTE MEASUREMENT BASICS (C40/NP80).....	36
SUBMICROLITER CELL MEASUREMENTS BASICS (C40).....	37
SAMPLE HANDLING TIPS .....	39
DATA TRANSFER .....	40
BATTERY OPERATION .....	41
<b>4. NANOPHOTOMETER APPLICATIONS.....</b>	<b>42</b>
NUCLEIC ACIDS .....	42
METHOD OVERVIEW.....	42

**MEASUREMENT PROTOCOL**.....42  
**CALCULATIONS** .....44  
**PROTEIN UV** .....48  
**METHOD OVERVIEW**.....48  
**MEASUREMENT PROTOCOL**.....49  
**CALCULATIONS** .....51  
**PROTEIN ASSAYS** .....55  
**METHOD OVERVIEW**.....55  
**MEASUREMENT PROTOCOL**.....56  
**CALCULATIONS** .....57  
**KINETICS** .....57  
**METHOD OVERVIEW**.....57  
**MEASUREMENT PROTOCOL**.....58  
**CALCULATIONS** .....59  
**OD600** .....59  
**METHOD OVERVIEW**.....59  
**MEASUREMENT PROTOCOL**.....61  
**CALCULATIONS** .....62  
**MORE APPS** .....62  
**MORE APPS: WAVELENGTH** .....62  
**METHOD OVERVIEW**.....62  
**MEASUREMENT PROTOCOL**.....62  
**CALCULATIONS** .....64  
**MORE APPS: WAVESCAN** .....64  
**METHOD OVERVIEW**.....64  
**MEASUREMENT PROTOCOL**.....64  
**CALCULATIONS** .....66  
**MORE APPS: ABSORBANCE RATIO** .....66  
**METHOD OVERVIEW**.....66  
**MEASUREMENT PROTOCOL**.....66  
**CALCULATIONS** .....68  
**MORE APPS: CONCENTRATION** .....68  
**METHOD OVERVIEW**.....68  
**MEASUREMENTS PROTOCOL**.....68  
**CALCULATIONS** .....70  
**MORE APPS: STANDARD CURVE** .....70  
**METHOD OVERVIEW**.....70  
**MEASUREMENTS PROTOCOL**.....70  
**CALCULATIONS** .....72  
**CUSTOM APPS** .....72  
**STORED RESULTS** .....72  
**STORED METHODS** .....73

<b>5. PREFERENCES</b> .....	<b>73</b>
GENERAL .....	73
<i>DATE AND TIME</i> .....	73
<i>DISPLAY</i> .....	74
<i>ABOUT</i> .....	74
<i>WiFi ACCESS POINT</i> .....	74
<i>ILLUMINATION SAMPLE WINDOW</i> .....	74
<i>ENABLE NANOVOLUME</i> .....	74
DYES 74	
WARNING MESSAGES .....	75
<i>BLANK CONTROL</i> .....	75
<i>SAMPLE QUALITY CONTROL</i> .....	76
NETWORK .....	76
<b>6. TROUBLESHOOTING</b> .....	<b>77</b>
SELF-CALIBRATION TEST .....	77
<b>7. ASSISTANCE</b> .....	<b>77</b>
SUPPORT .....	77
REPORT PROBLEM .....	77
SOFTWARE MAINTENANCE .....	78
<i>RESET</i> .....	78
<i>SOFTWARE UPDATE</i> .....	78
<i>CREATE LOG FILE</i> .....	79
<i>INSTALL PRINTER DRIVER</i> .....	79
LEAGAL .....	79
CONTACT IMPLEN .....	79
<b>8. MAINTENANCE</b> .....	<b>80</b>
MAINTENANCE FREE TECHNOLOGY .....	80
REPLACEMENT PARTS .....	81
CLEANING AND GENERAL CARE .....	81
MESSAGES .....	81
<i>IMPORTANT WARNING MESSAGES:</i> .....	82
<i>IMPORTANT ALERT MESSAGES:</i> .....	83
<b>9. WARRANTY</b> .....	<b>85</b>
<b>10. APPENDIX 1- INSTALACE SPEKTROFOTOMETRU</b> .....	<b>86</b>
BEZPEČNOSTNÍ INFORMACE .....	86
VYBALENÍ A UMÍSTĚNÍ .....	87
<b>11. APPENDIX 2 - INSTALLATIE SPECTROFOTOMETER</b> .....	<b>88</b>
VEILIGHEIDSINFORMATIE .....	88
UITPAKKEN EN PLAATSEN .....	89
<b>12. APPENDIX 3 - INSTALLATION DU SPECTROPHOTOMÈTRE</b> .....	<b>91</b>
INFORMATIONS DE SÉCURITÉ .....	91

DÉBALLAGE ET POSITIONNEMENT .....	92
<b>13. APPENDIX 4 – INSTALLATION SPECTROPHOTOMETER.....</b>	<b>94</b>
SICHERHEITSINFORMATIONEN.....	94
AUSPACKEN UND PLATZIEREN.....	94
<b>14. APPENDIX 5 – INSTALLAZIONE DELLO SPETTROFOTOMETRO .....</b>	<b>96</b>
INFORMAZIONI DI SICUREZZA.....	96
DISIMBALLAGGIO E COLLOCAZIONE.....	97
<b>15. APPENDIX 6 – INSTALACJA SPEKTROFOTOMETRU.....</b>	<b>98</b>
INFORMACJE DOTYCZĄCE BEZPIECZEŃSTWA.....	98
ROZPAKOWYWANIE I UMIEJSCOWIENIE .....	99
<b>16. APPENDIX 7 – INSTALACIÓN DEL ESPECTROFOTÓMETRO.....</b>	<b>101</b>
INFORMACIÓN DE SEGURIDAD .....	101
DESEMBALAJE E INSTALACIÓN .....	102
<b>17. APPENDIX 8 – SPEKTROFOTOMETERINSTALLATION .....</b>	<b>103</b>
SÄKERHETSINFORMATION .....	103
UPPACKNING OCH PLACERING .....	104
<b>18. APPENDIX 9 – INSTALAÇÃO DE ESPECTROFOTÓMETRO .....</b>	<b>105</b>
INFORMAÇÕES DE SEGURANÇA .....	105
DESEMBALAMENTO E POSICIONAMENTO .....	106
<b>19. ALPHABETICAL APPENDIX.....</b>	<b>108</b>

# 1. NANOPHOTOMETER AT A GLANCE

## NANOPHOTOMETER OVERVIEW

The Implen NanoPhotometer spectrophotometer is a mobile, simple to use UV/Visible instrument with a CCD array detector with options ranging from nanovolume to standard cuvette and all-in-one solutions. The N50, N60, and NP80 spectrophotometers measure sample volumes ranging from 0.3 - 2  $\mu$ l. The NP80 model also has the capability to measure samples using a standard cuvette. The C40 is a mobile standard cuvette solution with the option to upgrade to accommodate nanovolume measurements.

The NanoPhotometer runs on a Linux-based operating system (NPOS) that is designed for the use of pre-programmed and custom applications with a high degree of flexibility and processing power.

Sample Compression Technology™ provides easy sample handling which is independent of surface tension. This technology squeezes the sample between two quartz surfaces allowing for unmatched precision and accuracy without the need for dilutions. Combined with our True Path Technology™ the system offers lifetime accuracy and precision without the need for maintenance or recalibration.

**Note:** It is recommended to use a properly calibrated pipette with high-quality tips to ensure delivery of appropriate sample volumes for nanovolume sample applications.

The Sample Control™ is a quality control technology to identify air bubbles, sample impurities, turbidity, lint residues and potential contaminations. Sample Control™ monitors handling characteristics and sample quality in real time to ensure that the measured concentrations are reproducible and most precise.

The Blank Control™ gives a warning message for blanks with high background. The high background can be caused by the blank, blank buffer, or by residues from previous users which are the main cause for inaccurate readings. The Blank Control™ will protect the user from wasting time and precious sample on inaccurate readings caused by high background blanks or inappropriate cleaning. For more detailed information see also Technical Note #1 Blank Control™.

## NANOPHOTOMETER NP80

Item Number	Touchscreen	Battery Pack
NP80	-	-
NP80-Touch	+	-
NP80-Mobile	+	+

## NANOPHOTOMETER N60

Item Number	Touchscreen	Battery Pack
N60	-	-
N60-Touch	+	-
N60-Mobile	+	+

## NANOPHOTOMETER N50


Item Number	Touchscreen	Battery Pack
N50	-	-
N50-Touch	+	-

## NANOPHOTOMETER C40

Item Number	Touchscreen	Battery Pack
C40	-	-
C40-Touch	+	-
C40-Mobile	+	+

## INSTRUMENT REAR PANEL (NP80/N60/N50/C40)

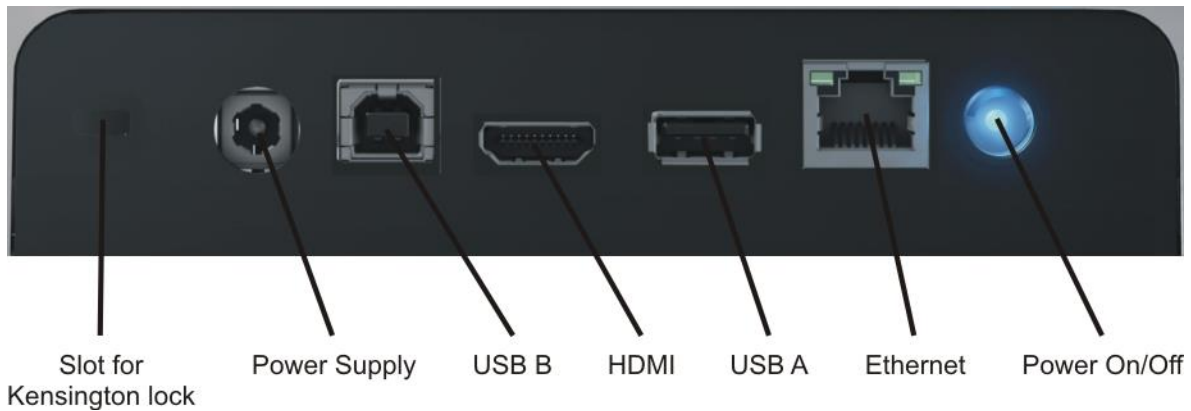## INSTRUMENT BOTTOM VIEW (NP80/N60/N50/C40)

Model name, device serial number and FCC ID are located on the identification plate on the bottom of the instrument.

## ACCESSORIES

### STANDARD ACCESSORIES

- **Connecting Cable**

USB connecting cable to connect the NanoPhotometer to a computer to control the NanoPhotometer via computer (please see page 21 Software Installation).

- **NanoPhotometer Power Adapter**

Power adapter for the NanoPhotometer and for charging the optional built-in battery pack.

**Note:** Use only the power adapter, supplied with your instrument or a replacement part from the manufacturer or your supplier.

- **Dust Cover**

### OPTIONAL ACCESSORIES

- **Didymium Glass Filter (C40/NP80)**

The certified didymium glass filter can be used to control the wavelength and photometric accuracy for the NanoPhotometer cuvette applications.

- **Standard Solution (NP80, N60, N50)**

The NanoPhotometer is recalibration free therefore it is not necessary to control the photometric accuracy on a regular basis. If the SOPs of a laboratory require a routine control of the photometric accuracy the standard solution can be used.

**Note:** Specifications of the standard solution are at least guaranteed for one year. Please see expiration date. Once the vial is opened it can be used for 30 minutes.

**Note:** Please read the Material Safety Data Sheet carefully prior to using this product.

- **IQ/OQ documentation (NP80, N60, C40)**

The NanoPhotometer IQ/OQ package consists of a Standard Solution and/or a Didymium Glass Filter and a corresponding Excel-based software tool depending on the type of instrument. The Excel-based software produces an automated report displaying all relevant data for an audit.

**Note:** The IQ/OQ documentation is not available for the NanoPhotometer N50.

- **Submicroliter Cell (C40 only)**

The submicroliter cell provides the option of upgrading the NanoPhotometer C40 to a device that handles nanovolume samples as well as cuvette samples. The sample volume range of the submicroliter cell ranges from a minimum of 0.3  $\mu\text{l}$  to 5.0  $\mu\text{l}$  with the path length options including 0.04 mm, 0.1 mm, 0.2 mm, 1 mm and 2 mm which correspond to virtual dilution factors of 1:250, 1:100, 1:50, 1:10, and 1:5 respectively.

- **DiluCell (C40/NP80 only)**

The DiluCell is a disposable cuvette with shortened path length for virtual dilution of cuvette-based samples. Due to the reduced path length the DiluCell provides an automatic dilution without the need of a physical dilution of higher concentration samples. There are two different sizes of DiluCells available DC 10 and DC 20 which allow an automatic 1/10 and 1/20 dilution of the sample respectively. Bypassing manual sample dilutions reduces dilution errors and cross-contamination making DiluCell ideal for GLP. Combined with small sample volume requirements and bubble free filling, the DiluCell allows for convenient spectrophotometric analysis from 340 - 950 nm.

- **Field Kit**

Implen aluminum roller case for safe transportation, designed to fit in most airplane overhead bins; special compartments for all necessary accessories, cleaning tools and samples.

**Note:** Delivered w/o NanoPhotometer and accessories


**Note:** Operate the NanoPhotometer only in the roller case when cover is open. Ensure that air circulation is possible. Switch the NanoPhotometer always off for transportation.

- **Dymo Label Printer**


It is possible to connect a Dymo Label printer to the NanoPhotometer for direct printing on self-adhesive labels. Recommended and tested printers are the Dymo LabelWriter 4XL (label size 10.3 x 15.8 cm) and the Dymo LabelWriter 450 (5.4 x 10 cm).

**Note:** The printouts are optimized for the label size of the Dymo LabelWriter 4XL. All other Dymo printers can be used. However, the font size will be zoomed to used paper size.

**Note:** Plug in the Dymo printer after starting the NanoPhotometer and the home screen is shown. Wait at least for 30 sec. for driver installation.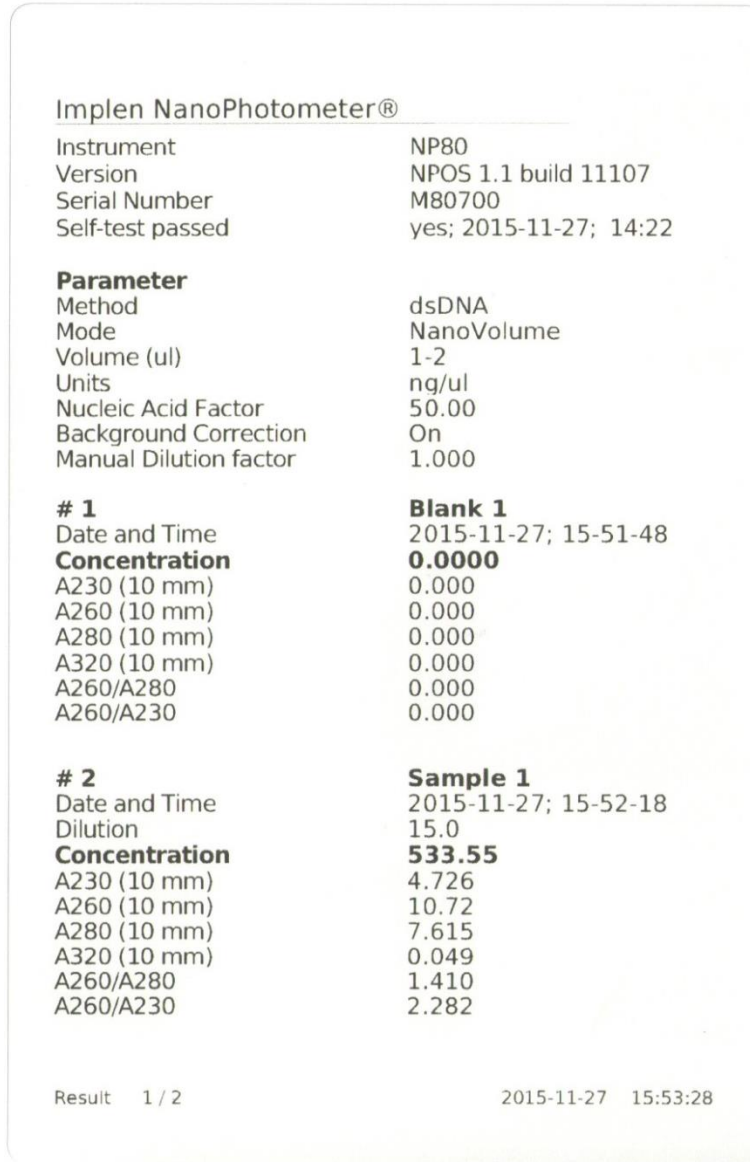

Printout for Dymo LabelWriter 4XL

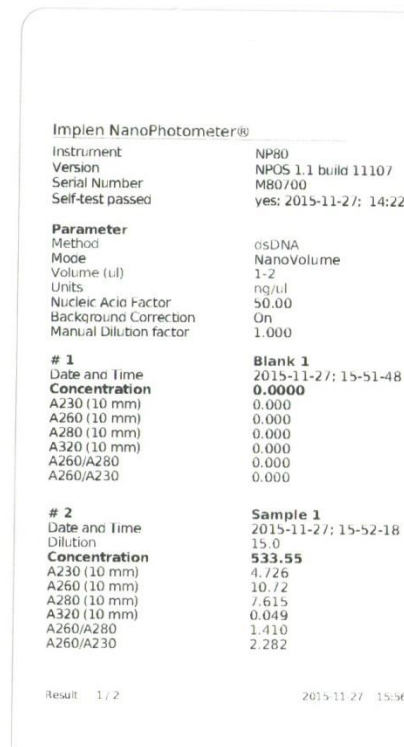Printout for Dymo LabelWriter 450

▪ **HP Printer**

It is possible to connect HP printer to the NanoPhotometer via USB cable for direct printout. The following HP printers are tested and released

- HP LaserJet 3030
- HP LaserJet m1522nf MFP
- HP LaserJet 400 color M451nw
- HP DeskJet 2543
- HP DeskJet 1110

Further HP printers are available upon request.

**Note:** Plug in the HP printer after starting the NanoPhotometer and the home screen is shown. Wait at least for 30 sec. for driver installation.

## CONNECTIVITY

### ▪ WiFi

The NanoPhotometer has a WiFi hotspot capability. This provides the option to control the NanoPhotometer by other WiFi devices like Windows computer, smartphones or tablets.

**SSID:** NanoPhotometer serial number; **password:** Implenuser

**Note:** Due to the limitations of some handheld devices saving to a wireless device is limited to 40 measurements per dataset. Larger datasets can be saved on the NanoPhotometer itself.

### ▪ USB A

There is a USB A port on the front and rear panel of the NanoPhotometer which is compatible with standard portable USB 2.0 storage devices for direct data transfer in a variety of formats including Excel. It is also possible to connect a mouse, keyboard, Dymo printer or HP printer directly to the NanoPhotometer.

**Note:** We recommend using FAT/FAT32 formatted 2.0 USB flash drives. Encrypted USB flash drives are not compatible.

**Note:** Cordless Bluetooth mice are not supported. Use only wired mice.

**Note:** Connect mouse and keyboard before starting the NanoPhotometer.

### ▪ USB B

There is a USB B port located on the rear panel of the instrument which is compatible with the USB cable provided to connect the NanoPhotometer to a computer. This USB connection can be used to control the NanoPhotometer via computer.

### ▪ LAN

There is an Ethernet (LAN) connection port on the rear panel of the instrument which enables the NanoPhotometer to connect with computers within close proximity. This Ethernet connection can be used to transfer data from the NanoPhotometer to a computer or to control the NanoPhotometer via a Windows computer. To access the directory of the NanoPhotometer from a Windows computer via LAN enter the serial number or IP (\\serial number / \\IP) of the NanoPhotometer in the Windows-Explorer and confirm with the enter key. For Mac OS X open the "Connect to Server" dialog in the "Go" menu of the MAC OS X Finder. Enter the instrument serial number or the active NanoPhotometer IP address in the server address field to connect.

**Note:** Plug in the LAN cable before starting the NanoPhotometer.

**Note:** The maximal LAN cable length is 10 meters. Bit rate is 1 Gbit/s.

### ▪ HDMI

There is an HDMI port located on the rear panel of the NanoPhotometer which is compatible with HDMI 1.4 cables (or better) to connect the NanoPhotometer to HDMI compatible monitors.

**Note:** The maximal HDMI cable length is 5 meters.

## NANOPHOTOMETER SPECIFICATIONS

### NanoVolume Performance

Detection Range dsDNA	1 ng/μl to 16,500 ng/μl (N50: 5 ng/μl to 7,500 ng/μl)
Detection Range BSA	0.03 mg/ml to 478 mg/ml (N50: 0.15 mg/ml to 217 mg/ml)
Minimum Sample Size	0.3 μl
Photometric Range (10 mm equivalent)	0.02 - 330 A (N50: 0.1 – 150A)
Path length	0.67 and 0.07 mm
Dilution Factor	15 and 140
Vortexer	2,800 rpm; tube size up to 2.0 ml

### Cuvette Performance

Detection Range dsDNA	0.1 ng / μl to 130 ng / μl
Detection Range BSA	0.003 mg / ml to 3.7 mg / ml
Photometric Range	0 - 2.6 A
Center Height (Z-Height)	8.5 mm
Cell Types	outside dimension 12.5 x 12.5 mm
Heating	37°C ± 0.5°C

### Optical Specifications

Wavelength Scan Range	200 – 900 nm (N50: 200 - 650 nm)
Time for Full Scan Range	3.5 – 6.0 seconds
Wavelength Reproducibility	± 0.2 nm (N50: ± 1nm)
Wavelength Accuracy	± 0.75 nm (N50: ± 1.5 nm)
Bandwidth	1.8 nm (N50: 5 nm)
Stray Light	< 0.5% at 240 nm using NaI (N50: < 2%) and < 1% at 280 nm using Acetone (N50: < 2%)
Absorbance Reproducibility	< 0.002 A (0.67 mm path) @ 280 nm (N50: < 0.004 A (0.67 mm path) @ 280 nm)
Absorbance Accuracy	< 1.75 % @ 0.7 A (0.67 mm path) @ 280 nm of the reading
Zero Stability	±0.003 A/hour after 20 min warm up @ 280 nm
Noise	0.002 A rms at 0 A @ 280 nm; 0.002 A (pk to pk) at 0 A @ 280 nm
Optical Arrangement	1 x 3648 CCD Array (N50: 1 x 1024 CCD Array)
Lamp	Xenon flash lamp
Lifetime	10 <sup>9</sup> flashes, up to 10 years

### Processing Power and Compatibility

Operating System	Linux based OS
Onboard Processor	Quad Core 1 GHz
Internal Storage	8 GB
Control Options	Onboard with built-in touchscreen, computer, smartphone, and tablet
Software Compatibility	Windows 7 (32 and 64 bit), Windows 8 (32 and 64 bit), OS X, iOS and Android
Min. Requirement	4 inch screen;
Smartphone/Tablet	Apple: iPad 2, iPhone 5 and iOS 6 Android version 4.4, Quad core 1.2 GHz with 1 GB RAM

### General Specifications

Size	200 mm x 200 mm x 140 mm
Weight	3.8 – 5.2 kg depending on configuration
Operating Voltage	90-250 V, 50/60 Hz, 60 W (90 W with battery pack), 18/19 VDC

Display	1024 x 600 pixels; touchscreen glove compatible
Certifications	CE, IEC/EN 61010-1:2012 and EN 61326-1:2013
Battery Pack	optional rechargeable lithium ion battery; 95Wh, 6,6Ah; Operation time: up to 8 h; min. charging cycles: 800
Battery Certification	IEC 62133 and UN38.3 transport test
In- and Output Ports	2x USB A, USB B, HDMI, Ethernet, WLAN
Additional Data Input	Mouse and keyboard options
Security	Slot for Kensington lock

Features and specifications are subject to change without notice.

US Patents 20080204755 and 20080106742

Windows is a trademark of Microsoft. Mac OS & iOS are trademarks of Apple. Android OS is a trademark of Google. Linux is a trademark of Linus Torvalds.

## 2. GETTING STARTED

### SPECTROPHOTOMETER INSTALLATION

**Note:** Translations of the Safety Information and Unpacking and Positioning are available in Appendix 1 (Czech), Appendix 2 (Dutch), Appendix 3 (French), Appendix 4 (German), Appendix 5 (Italian), Appendix 6 (Polish), Appendix 7 (Spanish), Appendix 8 (Swedish) and Appendix 9 (Portuguese)

#### SAFETY INFORMATION

Before commencing installation, please take time to familiarize yourself with warning labels and symbols on your instrument and their meaning. These are to inform you where potential danger exists or particular caution is required. Improper use may cause personal injuries or damage to the instrument. The instrument must only be operated by appropriately trained and experienced personnel. Please read the complete user manual prior to use.

 direct current

Do not open the instrument as this can expose the operator to electrical power, UV light, and delicate fiber optics or damage the instrument.

If the submicroliter cell is used with the NanoPhotometer C40 make sure that the lid is placed on the measurement head prior to the measurement.

**WARNING:** UV exposure when measuring without lid.

Do not use damaged power cords, accessories, and other peripherals with your NanoPhotometer. Use only the delivered and specified power supply/charger.

Do not expose the NanoPhotometer to strong magnetic, electrical fields, water, chemicals or any type of liquid as heavy rain or moisture.

Do not put the instrument into fire, as it may swell or explode (battery). Do not store at or use near any type of heat source, especially temperatures above 60°C or in an explosive atmosphere.

Do not leave your NanoPhotometer on your lap or near any part of your body to prevent discomfort or injury from heat exposure.

Do not place objects on top of the NanoPhotometer.

The NanoPhotometer with battery pack (mobile version) has to be switched off during transportation. The on/off button has to be protected during transportation from turning on itself caused by shock or vibration.

Biological samples may contain or have the potential to transmit infectious diseases. Be aware of the health hazard presented by such samples and wear appropriate protective equipment. Handle such samples with the greatest of care and according to applicable regulatory and organization requirements before working with such potential infectious materials.

**Note:** Do not spill any biological samples on instrument components. If spill occurs, disinfect the instrument immediately following your laboratory protocols and the cleaning instruction of the instrument (see page 80 Maintenance).

The symbol  on the product, or on the documents accompanying the product, indicates that this appliance may not be treated as household waste. Instead it shall be handed over to the applicable collection point for the recycling of electrical and electronic equipment. Disposal must be carried out in accordance with local environmental regulations for waste disposal.

## UNPACKING AND POSITIONING

**Note:** Translations of the Safety Information and Unpacking and Positioning are available in Appendix 1 (Czech), Appendix 2 (Dutch), Appendix 3 (French), Appendix 4 (German), Appendix 5 (Italian), Appendix 6 (Polish), Appendix 7 (Spanish), Appendix 8 (Swedish) and Appendix 9 (Portuguese)

Check the contents of the package against the delivery note. If any shortages are discovered, inform your supplier immediately.

Inspect the instrument for any signs of damage caused in transit. If any damage is discovered, inform your supplier immediately.

Ensure your proposed installation site conforms to the environmental conditions for safe operation: indoor use or dry environment.

**Note:** Do not expose your NanoPhotometer near liquids, chemicals, rain, moisture or dusty environments.

Temperature range 10 - 40°C; If the cuvette heating is used the range is 10 - 30°C.

If the instrument is subjected to extreme temperature changes, it may be necessary to allow the instrument to equilibrate. Turn the instrument off and then on again once thermal equilibrium has been established (~2 - 3 hours).

Maximum relative humidity (non-condensing) of 80% and up to 31°C decreasing linearly to 50% at 40°C.

The instrument must be placed on a stable, level surface that can support 4 - 5 kg and so that air can circulate freely around the instrument. Ensure while powered on that no materials reduce air circulation. Avoid direct sunlight as it may bleach parts of the instrument and can cause damage to plastic parts.

The equipment should be positioned such that in the event of an emergency the main plug can be easily located and removed.

Carry the instrument always by holding the main corpus of the instrument and not e.g. on the optional attached display or NanoVolume pedestal.

The equipment must be connected to power with the 60W power supply/cord supplied by Implen (90W with battery pack). The power outlet must have a protective conductor (earth/ground). It can be used on 90 - 250 V, 50 - 60 Hz power supply system.

For Mobile models please load the battery pack at least for 3 hours prior to the first use.

Turn the instrument on using the power button on the rear panel after it has been plugged in. The instrument will perform a series of self-diagnostic checks.

Please contact original supplier immediately if technical or sample handling difficulties are experienced.

**Note:** If this equipment is used in a manner not specified or in environmental condition not suitable for safe operation, the protection provided by the equipment may be impaired and the instrument warranty voided.

## SOFTWARE INSTALLATION

### NPOS OVERVIEW

NPOS is a Linux-based operating system designed for the NanoPhotometer.

NPOS can store data either to a common directory or be configured to save to independent directories according to file format and/or instrument.

NPOS can save data in an Implen .IDS format, PDF or as an Excel format file.

**Note:** PDF and Excel files cannot be opened on the NanoPhotometer. The files need to be transferred to a computer or device where Excel or a PDF reader is installed.

**Note:** Please do not connect the instrument to a computer until the NanoPhotometer NPOS software is installed on the computer.

### REQUIREMENTS AND COMPATIBILITY

The NPOS user interface is designed that all features can be operated by using a touchscreen. If the software is installed on a computer without touchscreen, the user interface can be operated by using a keyboard and a mouse. Before starting the installation process ensure the device is compatible with the requirements for installation.

#### Compatible Control Devices

##### Computer:

PC: Windows 7 / Windows 8 (32 bit or 64 bit)

Mac: OS X

##### Tablets (minimum requirements):

iPad 2: iOS6

Android (quad core 1.2 GHz with 1 GB RAM): Android version 4.4

**Smartphones (minimum requirements):**

iPhone 5: iOS6

Android (quad core 1.2 GHz with 1 GB RAM): Android version 4.4

Windows is a trademark of Microsoft. Mac OS & iOS are trademarks of Apple. Android OS is a trademark of Google. Linux is a trademark of Linus Torvalds.

**Note:** There are two user interfaces of the software available; one for built-in touchscreen, computer and tablets and one for smartphones.

## INSTALLING SOFTWARE ON COMPUTER

The NanoPhotometer software can be installed on compatible PC and Mac computer systems. Various operating systems and PC hardware may cause the set-up procedure to differ from that described here. This process is given as guidance only; it may need adaptation for other systems.


**Note:** Do not connect the NanoPhotometer to the PC/MAC before NPOS installation.

**Note:** If the software is already installed on the PC, remove the USB cable and uninstall the NPOS software first.

**Note:** The Windows and Mac installation file is provided on the [Implen](#) USB flash drive which is included in the NanoPhotometer delivery and are available for free download in the download section of the [Implen](#) webpage.

### ▪ NPOS installation for single user on Windows computer

1. Launch the NPOS installation file (NanoPhotometer.msi) and follow the installation routine. The installer should be logged with full administration rights. If you have insufficient privileges, installation may fail. If in doubt consult your PC administrator.
2. Open the NPOS software and choose the preferred connection. For USB connect the NanoPhotometer with the provided USB cable to the PC, for WiFi connection ensure stable WiFi connection between PC and NanoPhotometer (**SSID:** Serial number, **password:** Implenuser) and for LAN connect the NanoPhotometer to the local network via Ethernet cable.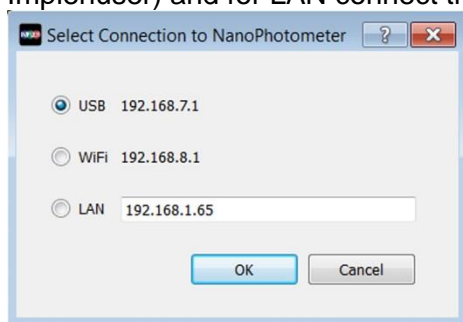


**Note:** If there is Avira installed on your computer, it is recommended to switch off the browser safety. This may interfere with the NPOS running on your computer.

### ▪ NPOS installation for multi user on Windows computer


1. Launch the NPOS installation file (NanoPhotometer.msi) and follow the installation routine.

Choose on the Licence Agreement dialog the option “Advanced” and on the following dialog “Install for all users of this machine”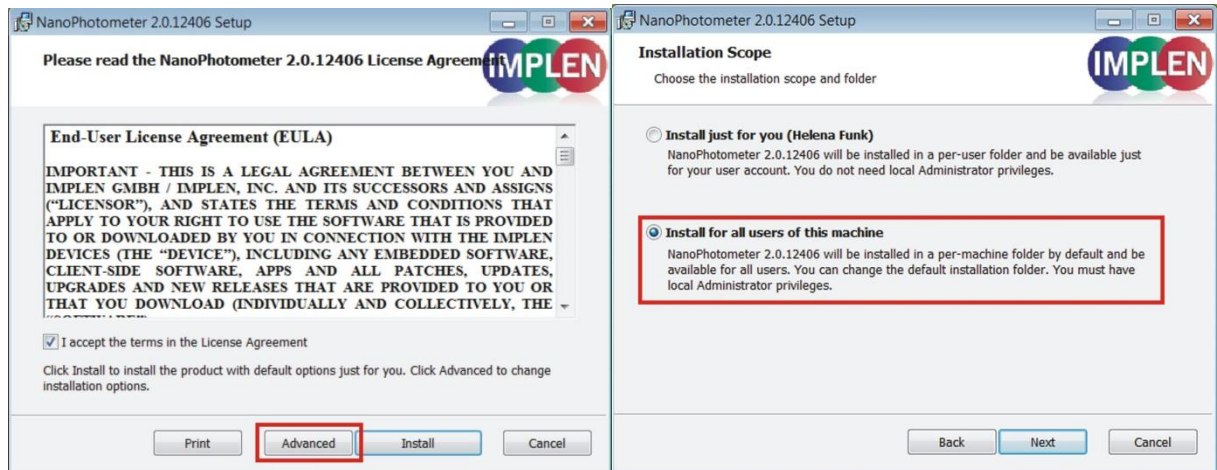

The installer should be logged with full administration rights. If you have insufficient privileges, installation may fail. If in doubt consult your PC administrator.

- Open the NPOS software and choose the preferred connection. For USB connect the NanoPhotometer with the provided USB cable to the PC, for WiFi connection ensure stable WiFi connection between PC and NanoPhotometer (**SSID**: Serial number, **password**: Implenuser) and for LAN connect the NanoPhotometer to the local network via Ethernet cable.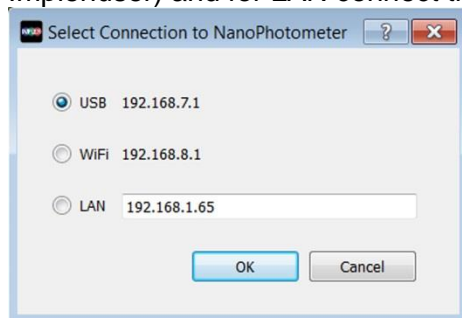

**Note:** If there is Avira installed on your computer, it is recommended to switch off the browser safety. This may interfere with the NPOS running on your computer.

## INSTALLING NANOPHOTOMETER APP ON TABLET OR SMARTPHONE

The NanoPhotometer App can be installed as an application on tablets and smartphones with compatible Android and iOS operating systems. The NanoPhotometer App is available for free download in the app store (Apple Store and Google Store).

- Download and install the NanoPhotometer App from the app store
- Connect the tablet or smartphone via WiFi to the NanoPhotometer.  
**SSID**: Serial number, **password**: Implenuser
- Open the NanoPhotometer App
- When connected via WiFi the NanoPhotometer will recognize the tablet as a remote control device and measurements can be initiated from the tablet or smartphone.
- Measurement results will be shown to the tablet or smartphone once taken

**Note:** In order to install the NanoPhotometer App on a tablet or smartphone, the device must have an established internet connection to access the app store for app download.

**Note:** The version of the app and the software of the NanoPhotometer should be the same. Different versions may have not full functionality.

## FIRST STEPS AND CONFIGURATION WIZARD

Starting the Implen NPOS the first time an Implen configuration wizard is shown. Please confirm the End User License Agreement (EULA) and select the country in which the NanoPhotometer is used and confirm.

## PRINTER INSTALLATION

For NanoPhotometer with software version 1.1.11147 or higher NO driver installation is necessary.

1. Switch NanoPhotometer on / home screen
2. Connect Dymo/HP printer via USB cable
3. Dymo/HP printer is ready to use after 30 seconds

For NanoPhotometer with previous software versions printer driver installation is necessary.

1. Switch NanoPhotometer on / home screen
2. Update to newest firmware version

### Update Process:

- a. Copy NPOS.bin to a USB flash drive into the root folder
- b. Insert the USB flash drive into your NanoPhotometer
- c. Select: "Assistance / Software Maintenance"
- d. Click on "Update"
- e. Wait until the NanoPhotometer reboots

**Note:** Update always first the firmware on the NanoPhotometer and in a second step the client software (PC/MAC, tablet, smartphone).

3. Go to "Assistance / Software Maintenance"
4. Start driver installation by a click to the "Install Driver for Printers" button
5. Wait until the following message is shown: "Installed driver for printer successfully."
6. Go to home screen
7. Connect Dymo/HP printer via USB cable
8. Wait for 30 seconds
9. Dymo/HP printer is ready to use

**Note:** Connect the printer after starting the NanoPhotometer or in the home screen. It is not possible to operate more than one printer at a time.


### 3. NANOPHOTOMETER BASICS

The NanoPhotometer product line offers a complete solution for Nanovolume (NP80/N60/N50/C40) and standard cuvette (C40/NP80) applications. With the NanoPhotometer NanoVolume applications the sample volume ranges from a minimum of 0.3  $\mu\text{l}$  to a maximum sample volume of 2.0  $\mu\text{l}$  (5  $\mu\text{l}$  for using the submicroliter cell/C40). Standard cuvette applications can be performed with 10 mm, 5mm, 2 mm and 1 mm path length quartz, glass, or plastic cuvettes with a center height of 8.5 mm.

#### APPLICATIONS OVERVIEW

The NanoPhotometer comes with pre-programmed applications as well as the ability to create custom applications. To select a method click/tap on the corresponding icon and the method will open immediately.

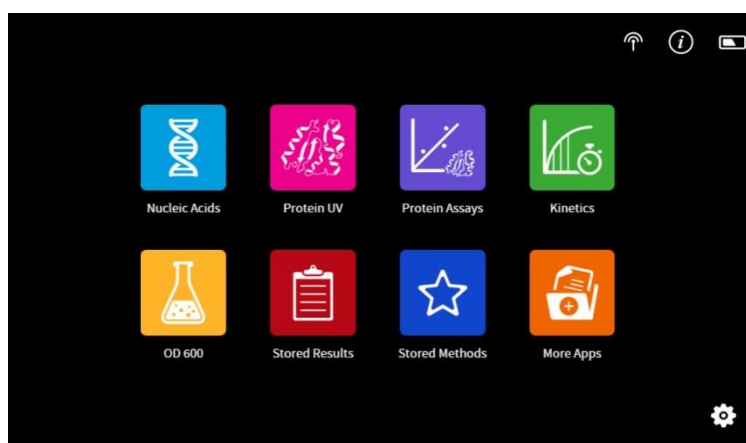#### Method Icons

**Nucleic Acids**

**Protein UV**

**Kinetics**

**Protein Assays**

**OD600**

#### Description

Concentration, purity, and dye incorporation for DNA, RNA, Oligo, and other nucleic acids

Protein UV determination at 280 nm (or in a range of 200 - 330 nm), purity and dye incorporation

Time vs. Absorbance readings

BCA (562 nm), Bradford (595 nm), Lowry (750 nm), and Biuret Assays (546 nm)

Measures cell density at 600 nm (or in a range of 200 – 900 nm)

- 
**Stored Results**    Archive of stored results
- 
**Stored Methods**    Collection of stored custom methods
- 
**More Apps**    Additional applications found on a secondary method screen
- 
**Wavelength**    Define one or multiple wavelength between 200 – 900 nm (N50: 200 – 650 nm) for absorbance measurements
- 
**Wavescan**    Define desired full scan range anywhere between 200 - 900 nm (N50: 200 - 650 nm)
- 
**Concentration**    Define extinction coefficient for automatic concentration calculations
- 
**Absorbance/ Ratio**    Define two wavelengths absorbance/ ratio calculation
- 
**Standard Curve**    Create a standard curve at a defined wavelength
- 
**Custom Apps**    Optional custom applications for personalized methods tailored to individual spectroscopy needs

## ICONS

Icon	Name	Action
	WiFi Hotspot	WiFi hotspot enabled; can be disabled in preferences
	Assistance	Opens the assistance page
	Battery Status	Shows the actual battery status (only shown with optional battery pack)
	Preferences	Opens the preferences page
	Home Screen	Returns to the home screen with application icons for method selection.
	Leave Method	Returns to the previous application selection
	Back	Returns to the previous page (smartphone only)
	Next/Confirm	Confirms parameter and opens the next screen (smartphone only)
	Delete	Deletes added functions in parameter; empties input windows
	Delete Data	Opens a delete dialog pop up
	Email Data	Opens an email dialog pop up (computer and tablet versions only)
	Save Data	Opens a save dialog pop up
	Print Data	Opens a print dialog pop up (only shown when Dymo/HP printer is successfully installed and connected)
	Add Folder	Adds a new folder to the directory
	Manage Data	Opens a dialog pop up with several action options including delete, rename or import folders/files/data as well as copying or move folders/files/data to defined directories
	Store Method	Opens a dialog pop up with the possibility to store the actual method parameter to a custom method
	Full Scale	Restores graph to original size without zoom
	Cancel	Returns to the previous screen without implementing any changes

## BUTTONS

### BLANK

Prior to any sample measurement at the start of a new method a blank measurement of either water or the buffer of the samples is required to give the NanoPhotometer a reference of what zero should be. It is recommended to re-apply the blank and measure it as a sample to ensure the graph of the blank spectra is a flat line.

### SAMPLE

To initiate the spectral scan of the sample press the sample button. The data will be temporarily stored until the method is exited at this time the user needs to define if the samples should be saved or deleted.

### Change to Cuvette

The Change to Cuvette button (NP80) is located at the bottom of the parameter window. The change to cuvette/nanovolume button toggles between nanovolume and cuvette applications. Selecting the Change to Cuvette button activates the cuvette compartment (opens the sliding cover and switches the red arrow LED on). Only cuvette measurements are possible when this mode is selected. Selecting the Change to Cuvette button while it is activated will deactivate the cuvette compartment (closes the sliding cover and switches the LED off) and only nanovolume measurements are possible.


**Note:** Please make sure that the nanovolume arm is closed during cuvette measurements.

### Change to NanoVolume

The Change to NanoVolume button (C40) is only shown when submicroliter cell applications are enabled in the preferences. The Change to NanoVolume/Cuvette button toggles between cuvette and submicroliter cell applications. With the NanoVolume submicroliter application the parameter of each method shows a dropdown list for lid selection (Lid 5, 10, 50, 100 and 250). Choose the correct lid and the results are automatically calculated considering the virtual dilution.

## SIDE TAB BAR

On the left side of the measurement screen there is a vertical tab bar that contains four tabs including: parameters, data, graph, and table. The different tabs allow organizing the measurement screen. It is possible to show or hide the different areas on the screen. Default screen for computer show all areas, for the built-in screen and tablet version the table is hidden.

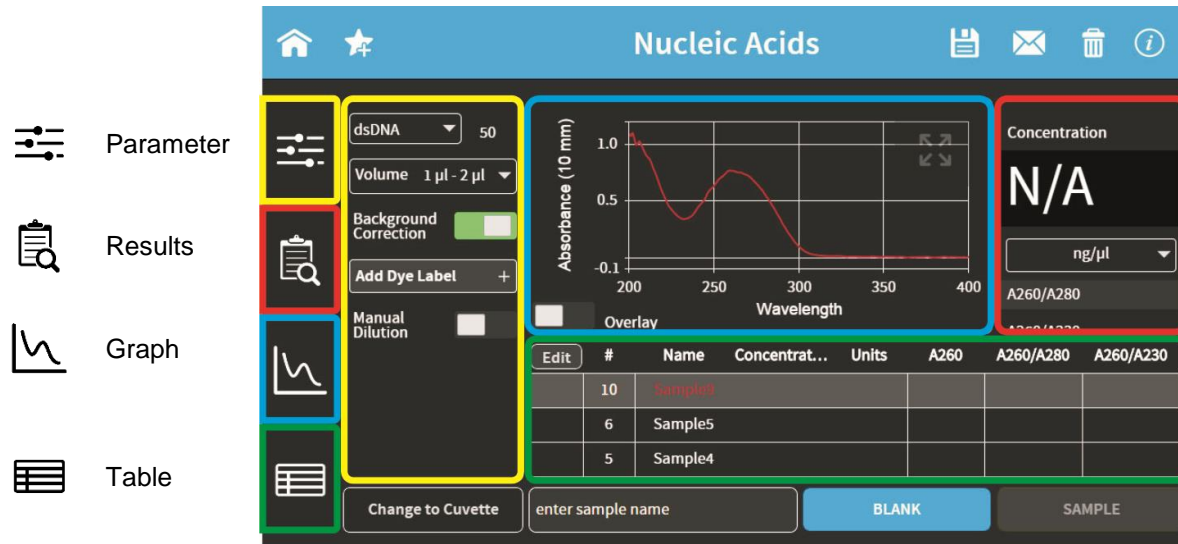**Note:** There is no tab bar available for smartphone versions. The parameters, results and graph screens are shown full screen. Parameters need to be confirmed (>) to get to the measurement screen. It is possible to toggle between the results and graph area by sliding left and right. There is no table area available on smartphones.

### Parameter area

In the parameter area it is possible to define all necessary parameters for a measurement as well as turn on cuvette mode and initiate cuvette heating. The default measurement screen includes the parameter area as open. It is automatically hidden when starting either a blank or sample measurement by clicking on the Blank or Sample button. It is also possible to hide the parameter area by tapping the parameter tab in the vertical side tab bar.

### Results area

The results area shows the method specific results of the latest measurement including concentration, absorbances, and relevant ratios from the last sample measured. It is also possible to change the units of the calculated concentrations in the results area with a drop-down selection menu.

**Note:** There are no data in the result view shown when the selection in the table is changed.

### ▪ Table area

The table area collects defined results for all samples completed during the active method. The first column of the table indicates whether the measurement is saved (📄) or not saved (blank field).

With the edit button the sample name of a single sample can be edited.

1. Select one sample in the table
2. Click on Edit Button
3. Change sample name
4. Confirm with the “Confirm” button

**Note:** The sample button is only active if one sample is selected in the table. Multi sample selection disables the Edit button.

### ▪ Graph area

The graph area shows a chart with the graph of the actual measurement or the selected line(s) in the table. There is an overlay toggle switch on the left bottom of the graph area. If the overlay option is enabled the graphs of the measurements will be automatically overlaid. To change the overlaid graphs select/deselect the samples in the table.

**Note:** It is only possible to overlay up to 30 graphs in a chart. If more than 30 data are selected a message will appear that says “More than 30 samples have been selected. Only 30 will be shown in graph.”

**Note:** The overlay button is not available on the NanoPhotometer and smartphones, only on tablets and computer versions.

It is possible to zoom in and out any area of the diagram (x- and y-axis). Undo the zoom by clicking the full scale icon (📏).

**Note:** Maximum zoom is 20 nm for the x-axis and 0.01A for the y-axis.

As legend option the sample name in the table is colored in the same color as the graph in the chart.

Clicking on the graph opens a pop up which shows the sample name, wavelength and absorbance of the selected wavelength. It is possible to show results of further graphs by changing the sample in the sample name drop-down.


Sample Name	L10
Wavelength	260
Absorbance	0.207

## DATA PROCESSING DIALOGS

### Print

Selecting the print data icon () opens a full screen overlay dialog window with various print options.

It is possible to connect Dymo or HP printer directly to the NanoPhotometer for printing via NanoPhotometer. On MAC and Windows computer it is also possible to use the printer connected to the computer or network for printing via Control Device.

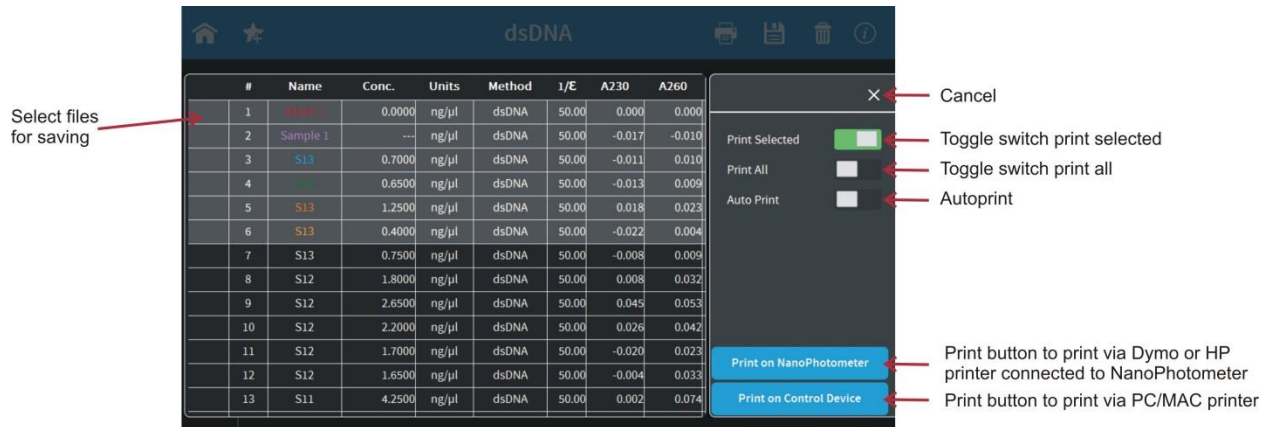**Note:** The print icon is shown, when a Dymo LabelWriter/HP printer is connected to the NanoPhotometer or in the computer version of the NPOS software.

**Note:** It is only possible to operate one printer at a time. Do not connect more than one printer to the NanoPhotometer.

#### Print selected

This option prints the current selection of data in the table. If the print selected toggle switch is green the print all toggle switch will automatically be deactivated.

#### Print all

The print all button (default setting) allows the user to print all of the data from the current method. If the print all toggle switch is green the save selected toggle switch will automatically be deactivated.

#### Auto print

If the auto print function is enabled, each measurement will be printed directly after the measurement on the connected printer.


**Note:** The default setting for auto print is off. If enabled in one method it is set to default on for all methods and needs to be switched off if not required.

**Note:** The auto print function is not available for printing via control device.

**Note:** The print option is not available in the smartphone apps.

## ▪ Save

Selecting the save data icon () opens a full screen overlay dialog window with various save options.

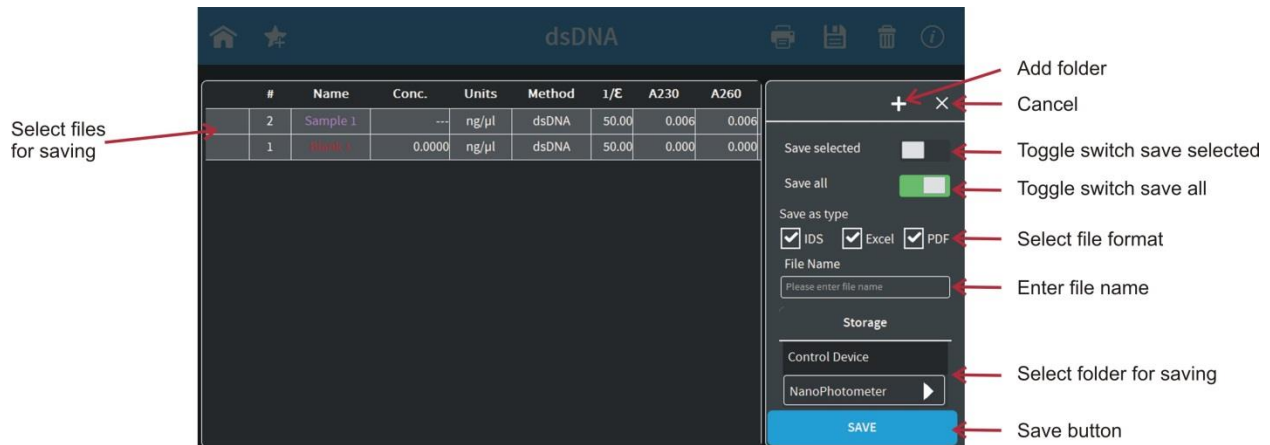### *Save selected*

This option saves the current selection of data in the table. If the save selected toggle switch is green the save all toggle switch will automatically be deactivated.

### *Save all*

The save all button (default setting) allows the user to save all data from the current method. If the save all toggle switch is green the save selected toggle switch will automatically be deactivated.

**Note:** On the smartphone app there is no save selected/save all toggle switch available; it is always the entire measurement session saved.

### *Save as type*

The save type option allows the user to specify in which file type the data are saved. The options of file types include Excel, PDF and Implen Document Source (IDS). It is possible to save different file formats in parallel.

**Note:** IDS files can only be saved on the NanoPhotometer and on a USB flash drive. It is not possible to save IDS files from opened data. In these cases the IDS tick box is greyed out.

**Note:** PDF and Excel files cannot be opened on the NanoPhotometer. The files need to be transferred to a computer or device where Excel or a PDF reader is installed.

**Note:** All data saved on the NanoPhotometer are saved on an internal micro SD card. We recommend to do data backups on a regularly basis to the hard drive of a computer or network. In the rarely case that the micro SD card crashes data loss cannot be ruled out.

## Implen Document Source File

Implen has a custom file format referred to as being an Implen Document Source (IDS) file. This is a hardcopy file to which no changes can be made. This file type contains all information

including raw data, results, values, and parameters. IDS files can only be opened with the Implen software.

### Excel File

Measurement data can be saved as an Excel file. This file format is compatible with Microsoft Excel software versions 2007/2010/2013 and can be edited. This file type contains all information including raw data, results, and parameters. The saved file contains only the selected information at the time the file was saved.

### PDF File

Measurement data can be saved as a PDF file. The saved file contains only the selected information at the time the file was saved.

#### *File Name*

Enter the name for the file. Allowed characters are: A...Z a...z 0...9 , . - ( ) @ ! = \_ ~ ; [ ] { } ‘

**Note:** Blank character is not allowed.

#### *Storage*

Selection option to choose the device the data should be saved to. The options include: NanoPhotometer, USB flash drive (if connected) and Control Device. If Control Device is selected the data will be transferred to the device that is currently connected with the NanoPhotometer.


**Note:** It is not possible to save IDS files to a computer (PC/Mac), tablets or smartphones.

### Autosave

To prevent data loss, all measurements are automatically stored as IDS file on the internal memory of the NanoPhotometer® right after each reading. These backup copies can be found in the Autosave folder on the device (StoredResults/NanoPhotometer/Autosave) for up to eight days. Files contain the base name Backup, the method name, and a time/date stamp.

- **Delete**

Selecting the delete icon (🗑️) opens a full screen overlay dialog window with delete options.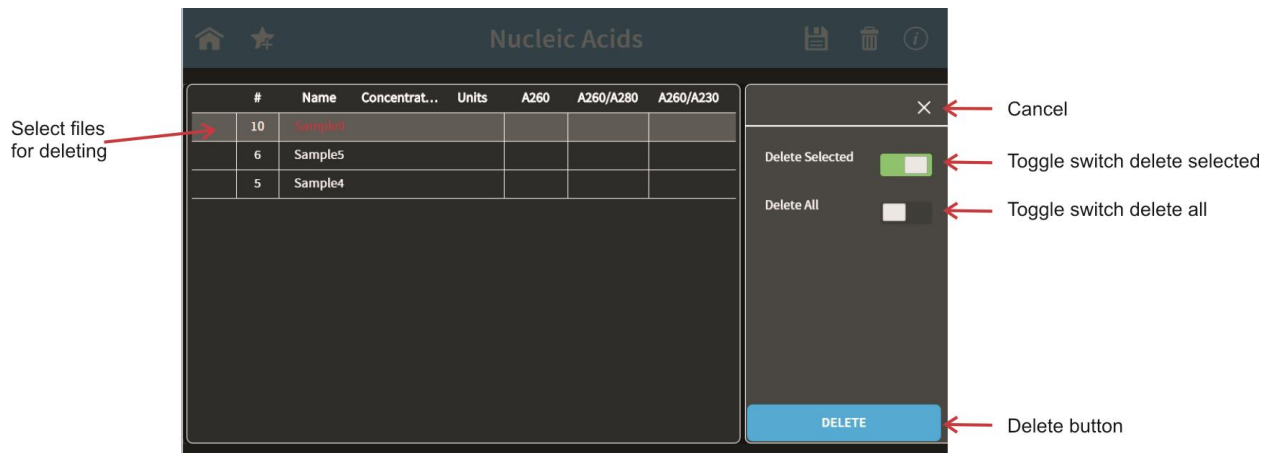

### Delete selected

This option deletes the current selection of data in the table. If the delete selected toggle switch is green the delete all toggle switch will automatically be deactivated.

### Delete all

The delete all button allows the user to delete all of the data from the current method. If the save all toggle switch is green the save selected toggle switch will automatically be deactivated.

Once the delete button is selected a message will appear asking “Do you want to delete all/selected files?” selecting cancel will return to the delete menu screen and selecting delete will delete the selected data.

**Note:** The delete function is not available for the software version designed for smartphones.


## BASIC OPERATION

The NanoPhotometer product line offers a complete solution for nanovolume (N50/N60/NP80/C40) and standard volume (C40/NP80) applications. With the NanoPhotometer nanovolume applications the sample volume ranges from 0.3  $\mu\text{l}$  to a maximum sample volume of 2.0  $\mu\text{l}$ . Standard Cuvette applications can be performed with 10 mm, 5 mm, 2 mm and 1 mm path length quartz, glass, or plastic cuvettes with a center height of 8.5 cm.

**Note:** NanoVolume application in the NanoPhotometer C40 is only available with the optional accessory “submicroliter cell” from Implen.

## NANOVOLUME MEASUREMENT BASICS (N50/N60/NP80)

1. Select a method depending on your sample and set the parameter for the measurement.
2. Ensure that the sample window and the mirror in the sample arm are clean.
3. Use the integrated vortexer (N60 and NP80 only) to mix your sample to achieve an homogenous sample

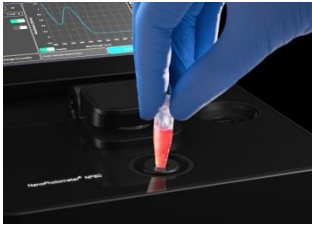4. Raise the sample arm and pipette the appropriate amount of blank solution onto the illuminated sample window. The illumination turns automatically off, lowering the arm.

**Note:** Do not overfill the well.  
**Note:** Save, low energy red light (LED)

5. Lower the sample arm and initiate a blank measurement with the blank button
6. Clean the measurement window and mirror on the sample arm with a slightly wet lint-free tissue. Use water, 70% ethanol or isopropanol.

**Note:** Make sure that the metal contact face (around the measurement window and the mirror) is clean.  
**Note:** Do not use aggressive solvents such as strong acids or bases or organic solvents at any time (see also page 40 solvent compatibility). If unsure please contact [support@implen.de](mailto:support@implen.de) for detailed information about your specific reagent/solvent.

7. It is possible to enter a sample name for each sample in the input window “enter sample name”.

**Note:** Allowed characters are: A...Z a...z 0...9 , . - ( ) @ ! = \_ ~ ; [ ] { } ‘  
**Note:** Blank character is not allowed.


8. Raise the sample arm and pipette the appropriate amount of sample solution onto the illuminated sample window. Upon completion of measurement raise sample arm, clean the surfaces and apply the next sample.  
**Note:** Parameter setting Volume 1 - 2  $\mu$ l adjusts the path length automatically. The

parameter setting Volume 0.3 µl measures only the 0.07 mm path length for higher concentrations (dsDNA > 420 ng/µl / BSA > 12.6 mg/ml).

**Note:** The sample window must be clean and residual fluff from any cleaning wipe must be removed for optimum performance.


## CUVETTE MEASUREMENT BASICS (C40/NP80)

The NanoPhotometer (NP80 and C40 only) is compatible with standard cuvettes having an 8.5 mm center height. The light path is indicated with a red status LED arrow for the NP80 model and with two white arrows for the C40 model.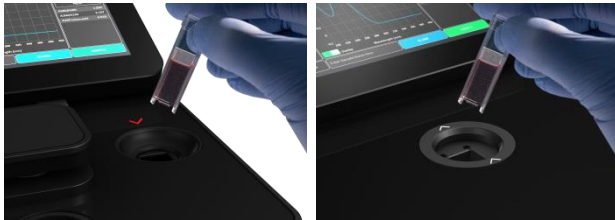

**Note:** For the NP80 the cuvette compartment needs to be activated by the “Change to Cuvette” button in the parameter area. Once the cuvette option is activated the sample compartment door will be opened automatically and a red arrow above the cuvette compartment will appear. The arrow indicates the light path.

1. Select a method depending on your sample and set the parameter for the measurement.
2. Open the cuvette cover by clicking the “Change to Cuvette” button below the parameter area (NP80).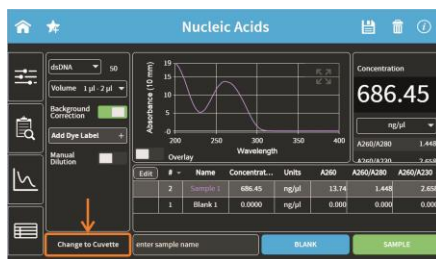

3. Add a blank solution to a cuvette and ensure that the filling volume is sufficient to allow light path to pass through the solution.
4. Insert cuvette into the cell holder.
5. Initiate a blank measurement with the blank button. Upon completion of measurement remove the cuvette.
6. Add sample to a cuvette and ensure that the sample volume is sufficient to allow light to pass through the sample.
7. It is possible to enter a sample name for each sample in the input window “enter sample name”.

**Note:** Allowed characters are: A...Z a...z 0...9 , . - ( ) @ ! = \_ ~ ; [ ] { } ‘  
**Note:** Blank character is not allowed.


8. Initiate a sample measurement with the sample button. Upon completion of measurement remove the cuvette.
9. Apply further samples.

## SUBMICROLITER CELL MEASUREMENTS BASICS (C40)

The Implen submicroliter cell can be used with the NanoPhotometer C40. Insert the submicroliter cell to the cell holder with the cell windows facing the light beam. We recommend facing the Implen logo to the right side. The light beam is directed from BACK to FRONT as indicated with the white arrows.

**Note:** For best performance insert the submicroliter cell always in the same direction.

1. Select a method depending on your sample.
2. Enable the NanoVolume application with the “Change to NanoVolume” button below the parameter area.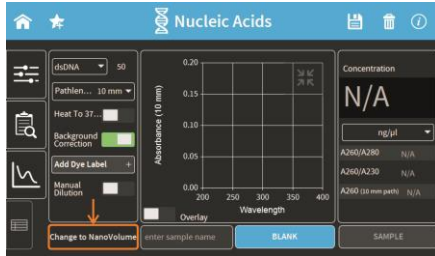

**Note:** If there is no “Change to NanoVolume” button available in the method go to preferences/general and enable the NanoVolume option. We guarantee our submicroliter cell specifications only for Implen submicroliter cells which are aligned to the NanoPhotometer C40 by Implen.

3. Set the parameter for the measurement. Choose the dilution lid depending on your sample concentration.

Lid	dsDNA (ng/µl)	BSA (mg/ml)	absorbance range 10 mm path
5 (optional)	2 – 375		0.05 – 7.5
10	5 – 750	0.15 – 22	0.10 – 15
50	25 – 3,750	0.70 – 108	0.50 – 150
100 (optional)	50 – 7,500	1.45 – 217	1.00 – 150
250 (optional)	125 – 18,750	3.63 – 543	2.50 – 375

**Note:** A new Blank is recommended when changing the dilution lid.

4. Ensure that the sample window and the mirror in the dilution lid are clean.
5. Pipette the appropriate sample volume (blank solution) onto the center of the measuring window. The required volumes are also shown in the parameter area of the software.

Lid	Sample volume	Path length	Dilution
5 (optional)	3.5 – 5 $\mu$ l	2 mm	1:5
10	1 – 3 $\mu$ l	1 mm	1:10
50	0.3 – 2 $\mu$ l	0.2 mm	1:50
100 (optional)	0.3 – 2 $\mu$ l	0.1 mm	1:100
250 (optional)	0.3 – 2 $\mu$ l	0.04 mm	1:250

**Note:** Do not overfill the well.

**Note:** It is a new blank for each change of the dilution lid recommended.

6. Ensure that for the measurements the lid fits exactly onto the positioning supports mounted to the body of the cell and initiate a blank measurement with the blank button.

**Note:** UV exposure. Do only initiate a measurement when the lid is closed.

7. Clean the measurement window and mirror in the lid with a slightly wet lint-free tissue. Use water, 70% EtOH or isopropanol.

**Note:** Do not use aggressive solvents such as strong acids or bases or organic solvents at any time (see also page 40 solvent compatibility). If unsure please contact [support@implen.de](mailto:support@implen.de) for detailed information about your specific reagent/solvent.

10. It is possible to enter a sample name for each sample in the input window “enter sample name”.

**Note:** Allowed characters are: A...Z a...z 0...9 , . - ( ) @ ! = \_ ~ ; [ ] { } ‘

**Note:** Blank character is not allowed.

8. Mix the sample very well to achieve a homogenous sample.
9. Apply the appropriate amount of sample solution onto the measurement window and initiate the sample measurement with the sample button. Upon completion of measurement remove the lid, clean the surfaces and apply the next sample.


## SAMPLE HANDLING TIPS

### ▪ Nanovolume Methods

- The NanoPhotometer includes an integrated vortexer (N60/NP80 only) to assure sample homogeneity. It is recommended to vortex every sample right before the measurement.
  - The sample window is illuminated (NP80/N60/N50 only) with a low energy red light to assist with accurate sample application. The red light is switched off once the sample arm is closed. It is possible to disable the illumination feature in preferences of the NPOS software.
  - The minimum volume that can be used for nanovolume samples is 0.3  $\mu\text{l}$  (starting for NP80/N60/N50 dsDNA > 420 ng/ $\mu\text{l}$  and BSA > 12.6 mg/ml and for submicroliter cell in the C40 dsDNA > 25 ng/ $\mu\text{l}$  and BSA > 0.7 mg/ml). For automatic path length setting at least 1  $\mu\text{l}$  is needed (NP80/N60/N50 only).
  - The maximum volume that can be used for nanovolume samples is 2.0  $\mu\text{l}$  (NP80/N60/N50) and 5  $\mu\text{l}$  for the submicroliter cell with lid 5.
  - The sample can be fully recovered after measurement with a pipette if desired.
- Note:** Minimal cross contamination cannot be avoided on a molecular level.
- Proper cleaning is important to ensure accurate measurements. In most cases a dry lint-free laboratory wipe is sufficient to clean the sample quartz surfaces. In the case of highly concentrated samples or certain proteins, the recommended procedure for cleaning is to use a slightly wet lint-free laboratory wipe (with water or 70% EtOH depending on sample type) to thoroughly clean the sample surface.
  - It is mandatory that the metal contact face around the measurement window and the mirror is clean.

### ▪ Cuvette Methods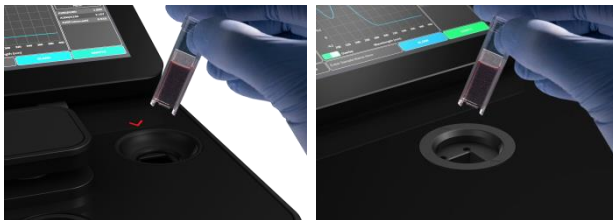

- The cuvette holder (NP80/C40 only) is compatible with standard 10 mm path length quartz, glass and plastic cuvettes with an optical height of 8.5 mm.
- It is also possible to use cuvettes with 5 mm, 2 mm or 1 mm path lengths, but there may be an adapter necessary. Please ask your cuvette supplier for suitable adapter.

- The minimum volume for accurate measurements depends on the cuvette type used; it is necessary that the light passes through the sample for accurate measurements. Center height is 8.5 mm.

**Note:** The cuvette holder is not removable. Do not pour any cleaning solution into the cuvette holder as larger amounts of liquids can get into the instrument and cause damage.

#### ▪ Solvent Compatibility (N50/N60/NP80/C40 with submicroliter cell)

Most solvents typically used in life science laboratories are compatible with the NanoPhotometer nanovolume sample surfaces. The following solvents are compatible for use with the NanoPhotometer models N50, N60, and NP80 at room temperature:

- Acetone ( $\leq 5\%$ )
- Acetonitrile
- Benzene
- Butanol
- Carbon tetrachloride
- Chloroform
- Ethanol
- Ether
- HEPES
- Hexane
- Isopropanol
- MES
- Methanol
- Methylene chloride
- MOPS
- Phenol ( $\leq 1\%$ )
- N-propanol
- Toluene
- Phosphate containing buffers
- PBS (pH 4-10)
- Citrate
- Borate
- Chloride salts
- Acids  $> \text{pH } 2$
- Bases  $< \text{pH } 10$

**Note:** Highly concentrated acids and bases are not recommended. It is recommended to wipe the sample surface with a lint-free laboratory wipe immediately upon completion of each measurement. For more information about compatibility of specific solvents not listed above, please contact the Implen support team ([support@implen.de](mailto:support@implen.de)) to check the compatibility.

## DATA TRANSFER

All data saved on the NanoPhotometer® can easily be accessed from and transferred to a computer within the lab through LAN, a USB cable or a password secured WiFi connection.

To access/transfer data from the NanoPhotometer Network Drive via LAN connection, the serial number or the NanoPhotometer IP needs to be entered in the address bar of the Windows Explorer (e.g. \\M80798\ or \\Assigned IP Address\). Using a USB cable connection \\192.168.7.1\ will provide access and for password secured WiFi connections \\192.168.8.1\ will need to be entered in the address bar to show the NanoPhotometer drive within Windows File Explorer. For Mac OS X open the "Connect to Server" dialog in the "Go" menu of the MAC OS X Finder. With a LAN connection enter the instrument serial number or the active NanoPhotometer IP address in the server address field to connect. For USB cable or WiFi connection use 192.168.7.1 or 192.168.8.1, respectively.

## BATTERY OPERATION

The optional battery operation is only available for the NanoPhotometer NP80/N60/C40 models (NP80-Mobile, N60-Mobile and C40-Mobile). A full battery pack can be operated for approximately 8 hours within normal usage, i.e. 20-30 measurements per hour. When the battery gets low there will be a first warning message “Battery low” and a beep is sounded. At this status it is possible to use the instrument under standard conditions at least for one hour. It is recommended to charge the battery at this status. If the battery is not charged there will be a second warning message when the battery is empty and several beep sounds before the instrument is automatically shutting down within 10 to 15 seconds.

**Note:** Unsaved data can be found in the auto save folder in Stored Methods. If the NanoPhotometer is shutting down during a measurement process the data of this measurement are lost.

The battery pack is designed for 800 full charging cycles. Afterwards the capacity and operation time of the battery may vary. A fully charged battery pack will self-discharge (switched off NanoPhotometer) over a time of 10 – 14 days. After this time recharging of the battery pack is necessary. Charging time for an empty battery pack is 3 hours.

## 4. NANOPHOTOMETER APPLICATIONS

The NanoPhotometer comes with pre-programmed applications as well as the ability to create custom applications. Any application method can be selected by tapping the icon once or clicking on the icon (computer based software).

**Note:** Selection of nanovolume or cuvette mode (NP80/C40) is within each method possible. The “Change to Cuvette” or “Change to NanoVolume” button is located at the bottom of the parameter area.

### NUCLEIC ACIDS

#### METHOD OVERVIEW


Nucleic acids in solution absorb light with a peak in the ultraviolet region of 260 nm. For determination of nucleic acid concentration in solution the absorbance at wavelength 260 nm is used along with the Beer-Lambert law. In addition to calculating concentrations of nucleic acids, absorbance measurements are also useful for estimating purity of nucleic acids by calculating the 260/280 nm and 260/230 nm ratios. Further, it is possible to determine the degree of labeling of nucleic acids with probes including fluorescent dyes.

The Sample Control™ gives useful information about air bubbles, sample impurities, turbidity, lint residues and potential contaminations. If the Sample Control™ detects any inconstancy an alert icon  is shown in the result/table area. A click on the alert icon shows additional information about the inconstancy.

#### MEASUREMENT PROTOCOL

1. Select the Nucleic Acids icon on the home screen.  
To change between NanoVolume and cuvette application (NP80/C40 only) use the Change to Cuvette/NanoVolume button below the parameter area.
2. Select the nucleic acid type in the drop-down of the parameter area.  
Options are: dsDNA, ssDNA, RNA miRNA, miRNA Sequence, Oligo, Oligo Sequence and Custom (see Table 1).
  - The miRNA sequence and Oligo sequence options allow having the sequence entered and the extinction coefficient will be automatically calculated.
  - Custom is set to 50 as a default and allows the extinction coefficient to be manually entered from the range of 15 - 150.

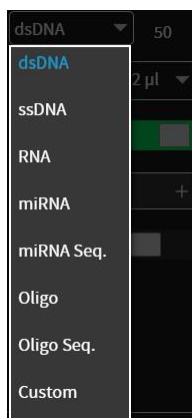3. **For NanoVolume application:**

Select the volume of sample to be applied.

**Note:** 1 - 2 µl (default): automatic path length change; 0.3 µl measures only the 0.07 mm path length (for samples with concentrations > 420 ng/µl dsDNA)

**For cuvette application:**


Select the path length depending on the used cuvette.

Options are: 1 mm, 2 mm, 5 mm and 10 mm

**For NanoVolume application with submicroliter cell (C40 only):**

Select Lid depending on your sample concentration for concentration ranges of different lids (see also page 37 Submicroliter Cell Measurements Basics). Required sample volume for each lid is shown below the lid selection dropdown.

**Note:** A new Blank is recommended when changing the dilution lid.

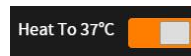
4. If it is desired to heat the sample to 37°C use the toggle switch to turn on cell holder heater. When the cuvette holder has 37°C the color changes to green.

**Note:** Only available for cuvette applications (NP80 and C40).

5. The background correction is enabled as the default. It is possible to disable the background correction with the toggle switch.

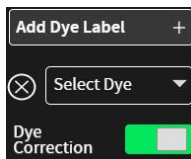6. If a dye labeled sample is used please add a dye by tapping on the Add Dye Label option and select a dye from the drop-down list. There is also an option for dye correction which can be en-/disabled with a toggle switch. Added Dyes can be deleted by tapping on the delete icon (⊗).

**Note:** If the used dye is not available in the drop-down list, please go to preferences and add a custom dye to the dye list.

7. Option to set/calculate a dilution factor for manual diluted samples.

8. Apply the blank ddH<sub>2</sub>O or buffer to the illuminated sample window for the reference measurement and select blank to initiate the reading.

**Note:** The illumination of the sample window can be switched off in the preferences.

9. Use a lint-free laboratory wipe to clean both the sample window and mirror prior to applying the next sample.

**Note:** It could be helpful to apply the blank a second time and read it as a sample to ensure a proper blank.

SAMPLE

- Apply sample to the sample window and press the sample button to initiate the measurement.

## CALCULATIONS

### ▪ Nucleic Acid Concentration

To determine the concentration of nucleic acids in solution, the absorbance as measured at the wavelength of 260 nm is used. The function describing the relationship between concentration and absorbance is a modification of the Beer-Lambert law equation. The concentration of the nucleic acid samples can be calculated with or without background correction depending on enabled/disabled background correction option.

#### Without background correction:

$$C = A_{260} * \epsilon_{nuc} * D$$

#### With background correction:

$$C = (A_{260} - A_{320}) * \epsilon_{nuc} * D$$

C	Concentration in ng/μl
A <sub>260</sub>	Absorbance at 260 nm (10 mm path)
A <sub>320</sub>	Absorbance at 320 nm (10 mm path)
D	Manual dilution factor
ε <sub>nuc</sub>	Extinction coefficient/nucleic acid factor in ng*cm/μl

**Table 1. Nucleic acids extinction coefficients (ε<sub>nuc</sub>)**

Type	ε <sub>nuc</sub>
dsDNA	50 ng*cm/μl
ssDNA	37 ng*cm/μl
RNA	40 ng*cm/μl
miRNA	33 ng*cm/μl
Oligo	33 ng*cm/μl
miRNA Seq.	calculated via extinction coefficient of constituent nucleotides entered
Oligo Seq.	calculated via extinction coefficient of constituent nucleotides entered
Custom	Option to enter any factor between 15 and 150 ng*cm/μl

### ▪ Dye-labeled Nucleic Acid Concentration

For dye-labeled nucleic acids, the concentration of the nucleic acid is calculated using a modified form of the Beer-Lambert equation. For these calculations, the instrument considers the absorption maximum of the dye and a certain dye-specific correction factor at 260 nm (see

Table 2). The concentration of a dye-labeled nucleic acid is calculated with or without background/dye correction as follows:

**With background and with dye correction:**

$$C = [(A_{260} - A_{320}) - (cf_{dye} * (A_{max,dye} - A_{320}))] * \epsilon_{nuc} * D$$

**With background and without dye correction:**

$$C = (A_{260} - A_{320}) * \epsilon_{nuc} * D$$

**Without background and with dye correction:**

$$C = [A_{260} - cf_{dye} * A_{max,dye}] * \epsilon_{nuc} * D$$

**Without background and without dye correction:**

$$C = A_{260} * \epsilon_{nuc} * D$$

C	Concentration in ng/μl
A <sub>260</sub>	Absorbance at 260 nm (10 mm path)
A <sub>320</sub>	Absorbance at 320 nm (10 mm path)
A <sub>max,dye</sub>	Absorbance value at the absorbance maximum of the dye (10 mm path)
ε <sub>nuc</sub>	Extinction coefficient/nucleic acid factor in ng*cm/μl
cf <sub>dye</sub>	Dye-dependent correction factor at 260 nm
D	Manual dilution factor

**Note:** Dye correction is only available for single dye selection.

▪ **Dye Concentration**

For dye-labeled nucleic acids, the concentration of the dye is calculated using a modified form of the Beer-Lambert equation. For these calculations, the instrument considers the absorption maximum of the dye, the dye-specific extinction coefficient (see Table 2). The dye concentration is calculated with or without background correction as follows:

**With background correction:**

$$C = \frac{(A_{max,dye} - A_{320}) * D}{\epsilon_{dye} * 10^6}$$

**Without background correction:**

$$C = \frac{A_{max,dye} * D}{\epsilon_{dye} * 10^6}$$

C	Concentration in ng/μl
---	------------------------

$A_{\text{max,dye}}$	Absorbance value at the absorbance maximum of the dye (10 mm path)
$A_{320}$	Absorbance at 320 nm (10 mm path)
$\epsilon_{\text{dye}}$	Extinction coefficient of dye in $\text{M}^{-1}\cdot\text{cm}^{-1}$
$D$	Manual dilution factor

### ▪ Frequency of Incorporation (FOI)

FOI is the degree of labeling based on dye incorporation in a labeled nucleic acid sample. It is generally expressed as the number of dye molecules incorporated per 1000 nucleotides. FOI can be calculated with or without background/dye correction as follows:

#### With background correction with dye correction:

$$\text{FOI} = \frac{324.5 * (A_{\text{max, dye}} - A_{320})}{\epsilon_{\text{dye}} * 10^6 * (A_{260} - A_{320} - \text{cf}_{\text{dye}} * (A_{\text{max,dye}} - A_{320}))}$$

#### With background correction without dye correction:

$$\text{FOI} = \frac{324.5 * (A_{\text{max, dye}} - A_{320})}{\epsilon_{\text{dye}} * 10^6 * (A_{260} - A_{320}) * \epsilon_{\text{nuc}}}$$

#### Without background correction and with dye correction:

$$\text{FOI} = \frac{324.5 * A_{\text{max, dye}}}{\epsilon_{\text{dye}} * 10^6 * (A_{260} - \text{cf}_{\text{dye}} * A_{\text{max,dye}}) * \epsilon_{\text{nuc}}}$$

#### Without background correction and without dye correction:

$$\text{FOI} = \frac{324.5 * A_{\text{max, dye}}}{\epsilon_{\text{dye}} * 10^6 * A_{260} * \epsilon_{\text{nuc}}}$$

FOI	Frequency of Incorporation (dye per 1,000 bases)
$A_{260}$	Absorbance at 260 nm (10 mm path)
$A_{320}$	Absorbance at 320 nm (10 mm path)
$A_{\text{max,dye}}$	Absorbance value at the absorbance maximum of the dye (10 mm path)
$\epsilon_{\text{dye}}$	Extinction coefficient of dye in $\text{M}^{-1}\cdot\text{cm}^{-1}$
$\epsilon_{\text{nuc}}$	Extinction coefficient/nucleic acid factor in $\text{ng}\cdot\text{cm}/\mu\text{l}$
$\text{cf}_{\text{dye}}$	Dye-dependent correction factor at 260 nm

**Table 2. Dye Types, Absorbance Max, Extinction coefficient, and dye-dependent correction factors**

NanoPhotometer Model	Dye Type	Absorbance maximum of Dye (nm)	Dye-dependent ext. coeff. $\epsilon_{\text{dye}}$ in $\text{M}^{-1}\cdot\text{cm}^{-1}$	Dye-dependent correction factor at 260 nm $\text{cf}_{\text{dye}}$
NP80/N60/N50/C40	Alexa Fluor 350	345	18,400	0.25
NP80/N60/N50/C40	Alexa Fluor 488	492	62,000	0.30
NP80/N60/N50/C40	Alexa Fluor 532	525	82,300	0.24
NP80/N60/N50/C40	Alexa Fluor 546	555	104,000	0.21
NP80/N60/N50/C40	Alexa Fluor 555	555	150,000	0.04
NP80/N60/N50/C40	Alexa Fluor 568	576	93,000	0.45
NP80/N60/N50/C40	Alexa Fluor 594	588	80,400	0.43
NP80/N60/C40	Alexa Fluor 647	650	239,000	0.00
NP80/N60/C40	Alexa Fluor 660	660	107,000	0.00
NP80/N60/C40	Alexa Fluor 680	680	164,000	0.00
NP80/N60/N50/C40	Cy3	550	150,000	0.08
NP80/N60/N50/C40	Cy3.5	581	150,000	0.08
NP80/N60/C40	Cy5	649	250,000	0.05
NP80/N60/C40	Cy5.5	675	250,000	0.05
NP80/N60/N50/C40	Oyster-500	503	78,000	0.29
NP80/N60/N50/C40	Oyster-550	553	150,000	0.05
NP80/N60/N50/C40	Oyster-556	560	155,000	0.03
NP80/N60/C40	Oyster-645	649	220,000	0.05
NP80/N60/C40	Oyster-650	653	200,000	0.04
NP80/N60/C40	Oyster-656	660	200,000	0.04
NP80/N60/N50/C40	Texas Red	593	85,000	0.23

#### ▪ Ratios

Reactions utilizing nucleic acids often require minimum purity standards. Common contaminants of nucleic acid samples include: proteins, organic compounds, and other. Based on the common contaminants of nucleic acid samples, the 260/280 and 260/230 ratios are calculated for nucleic acids to give an indication of the purity of the samples. Pure DNA and RNA preparations have expected 260/280 ratios of  $\geq 1.8$  and  $\geq 2.0$  respectively. An elevated absorbance at 230 nm can indicate the presence of impurities as well; 230 nm is near the absorbance maximum of peptide bonds and also indicates buffer contamination since TRIS, EDTA and other buffer salts absorb at 230nm. When measuring RNA samples, the 260/230 ratio should be  $> 2.0$ ; a ratio lower than

this is generally indicative of contamination with guanidinium thiocyanate, a reagent commonly used in RNA purification and which absorbs over the 230 - 260 nm range. If a ratio is detected out of the expeptable range an alert icon  is shown in the the results/table area. A click on the alert icon shows additional information.

The ratios are calculated with or without background correction according to if the background correction is activated during the measurements or not as follows:

#### Without background correction:

$$260/280 \text{ ratio} = \frac{A_{260}}{A_{280}}$$

$$260/230 \text{ ratio} = \frac{A_{260}}{A_{230}}$$

#### With background correction:

$$260/280 \text{ ratio} = \frac{A_{260} - A_{320}}{A_{280} - A_{320}}$$

$$260/230 \text{ ratio} = \frac{A_{260} - A_{320}}{A_{230} - A_{320}}$$

## PROTEIN UV

### METHOD OVERVIEW

The Protein UV method exploits the inherent absorbance of proteins at 280 nm in combination with the Beer-Lambert Law, where each protein is characterized by a protein specific extinction coefficient ( $\epsilon$ ) which can be used to determine total protein concentration of a solution. The intrinsic absorbance of proteins is due to the presence of aromatic amino acids in their structure, primarily tryptophan and tyrosine, as well as cysteine (oxidized cysteine residues in a disulphide bond). The aromatic amino acid residues in a protein containing tryptophan and tyrosine exhibit strong intrinsic absorbance at 280 nm, with a lesser contribution by phenylalanine. Therefore, it is the aromatic amino acid residues which dictate the extinction coefficient at 280 nm for a protein.

The most straightforward method to determine concentration of a purified, homogenous protein with a known extinction coefficient ( $\epsilon$ ) is by direct measurement of UV280 provided as long as the protein contains no prosthetic groups with strong absorption in the same region. However, for unknown proteins including homogenous protein mixtures, it is possible to make direct  $A_{280}$  measurements using a composite  $\epsilon$  value derived from comparison of many proteins, although this will only provide an approximate but close estimate of the true protein concentration.

The NanoPhotometer determines protein concentration by performing calculations based on specific  $\epsilon$  values, either pre-programmed in the instrument or entered manually by the user. Extinction coefficient ( $\epsilon$ ) values at 280 nm vary greatly for different proteins due to their

particular aromatic amino acid content. Fixed  $\epsilon$  values are pre-programmed in the software for certain proteins (see Table 3). However, if the protein of interest is not included in the pre-programmed methods it is also possible to manually enter the specific  $\epsilon$  for the protein of interest using the Mol. Ext. Coefficient or Ext. Coefficient protein option. For correct calculation, it is necessary to supply either: a) the mass extinction coefficient ( $\epsilon$  in  $l/g \cdot cm$ ); or b) the molar extinction coefficient ( $\epsilon_M$  in  $M^{-1} \cdot cm^{-1}$ ), and the molecular weight of the protein expressed in molar mass units (g/mol).

To determine the degree of dye labelling of a protein, the absorbance measured at the wavelength corresponding to the absorbance maximum of the fluorescence dye is used (see Table 3). The corresponding extinction coefficient of the dye is used along with the Beer-Lambert Law to determine the dye concentration.


**Note:** It is important to ensure the extinction coefficient and units entered are correct in order to ensure that calculations are performed properly for accurate concentration values.

The Sample Control™ gives useful information about air bubbles, sample impurities, turbidity, lint residues and potential contaminations. If the Sample Control™ detects any inconstancy an alert icon  is shown in the result/table area. A click on the alert icon shows additional information about the inconstancy.


## MEASUREMENT PROTOCOL

1. Select the Protein UV icon on the home screen.  
To change between NanoVolume and cuvette application (NP80/C40 only) use the Change to Cuvette/NanoVolume button below the parameter area.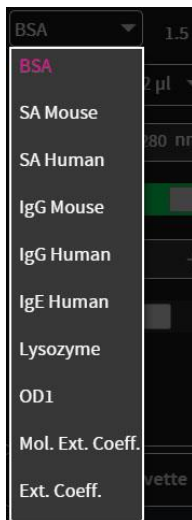

2. Select the protein type in the drop-down of the parameter area:  
Options are: BSA, SA Mouse, SA Human, IgG Mouse, IgG Human, IgE Human, Lysozyme, OD1, Mol. Ext. Coefficient, Ext. Coefficient.
  - For Mol. Ext. Coefficient enter Molecular Weight (default 66,400 g/mol) and Mol. Ext. Coefficient (default  $44,289 M^{-1} \cdot cm^{-1}$ )
  - For Ext. Coefficient enter Ext. Coefficient (default  $0.667 l/g \cdot cm$ )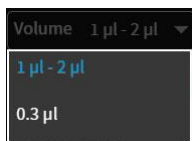

3. **For NanoVolume application:**  
Select the volume of sample to be applied.  
**Note:** 1 - 2  $\mu l$  (default): automatic path length change; 0.3  $\mu l$  measures only the 0.07 mm path length (possible for samples with concentrations e.g. BSA

> 12.6 mg/ml)

**For cuvette application:**

Select the path length depending on the used cuvette.

Options are: 1 mm, 2 mm, 5 mm and 10 mm

**For NanoVolume application with submicroliter cell (C40 only):**


Select Lid depending on your sample concentration for concentration ranges of different lids (see also page 37 Submicroliter Cell Measurements Basics).

Required sample volume for each lid is shown below the lid selection dropdown.

**Note:** A new Blank is recommended when changing the dilution lid.


- The background correction is enabled as the default. It is possible to disable the background correction with the toggle switch.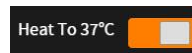

- If it is desired to heat the sample to 37°C use the toggle switch to turn on cell holder heater. When the cuvette holder has 37°C the color changes to green.

**Note:** Only available for cuvette applications (NP80 and C40).

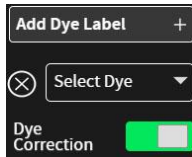- If a dye labeled sample is used please add a dye by tapping on the Add Dye Label option and select a dye from the drop-down list. There is also an option for dye correction which can be en-/disabled with a toggle switch. Added Dyes can be deleted by tapping on the delete icon (⊗).

**Note:** If the used dye is not available in the drop-down list, please go to preferences and add a custom dye to the dye list.

- Option to set/calculate a dilution factor for manual diluted samples.

- Apply the blank ddH<sub>2</sub>O or buffer to the illuminated sample window for the reference measurement and select blank to initiate the reading.

**Note:** The illumination of the sample window can be switched off in the preferences.

- Use a lint-free laboratory wipe to clean both the sample window and mirror prior to applying the next sample.

**Note:** It could be helpful to apply the blank a second time and read it as a sample to ensure a proper blank.

- Apply sample to the sample window and press the sample button to initiate the measurement.

## CALCULATIONS

### ▪ Protein UV280 Concentration

The protein concentration in the Protein UV method is calculated with the absorbance value of the sample at 280 nm along with the extinction coefficient defined by the user. The protein concentration is calculated with or without background correction as follows:

**With background correction:**

$$C = (A_{280} - A_{320}) * \epsilon_{\text{prot}} * D$$

**Without background correction:**

$$C = A_{280} * \epsilon_{\text{prot}} * D$$

C	Concentration (mg/ml)
A <sub>280</sub>	Absorbance at 280 nm (10 mm path)
A <sub>320</sub>	Absorbance at 320 nm (10 mm path)
ε <sub>prot</sub>	Extinction coefficient/protein factor in g*cm/l
D	Dilution factor

**Table 3. Protein extinction coefficients (ε<sub>prot</sub>)**

Type	ε <sub>prot</sub> [g*cm/l]	Ext. Coeff. [l/g*cm]	Mol. Ext. Coeff. [M <sup>-1</sup> *cm <sup>-1</sup> ]	MW [g/mol]
BSA	1.50	0.667	44,289	66,400
SA Mouse	1.49	0.670	44,220	66,000
SA Human	1.72	0.582	40,370	69,365
IgG Mouse	0.71	1.400	224,000	160,000
IgG Human	0.74	1.360	204,000	150,000
IgE Human	0.65	1.530	290,700	190,000
Lysozyme	0.38	2.640	37,984	37,984
OD1	1.00	N/A	N/A	N/A

### ▪ Dye-labeled Protein UV280 Concentration

For dye-labeled proteins, the concentration of the protein is calculated using a modified form of the Beer-Lambert equation. For these calculations, the instrument considers the absorption maximum of the dye, and a certain dye-specific correction factor at 280 nm (see Table 4). The dye concentration is calculated with or without background/dye correction for are as follows:

**With background and with dye correction:**

$$C = [A_{280} - A_{320} - cf_{\text{dye}} * (A_{\text{max,dye}} - A_{320})] * \epsilon_{\text{prot}} * D$$

**With background and without dye correction:**

$$C = (A_{280} - A_{320}) * \epsilon_{\text{prot}} * D$$

**Without background and with dye correction:**

$$C = (A_{280} - cf_{dye} * A_{max,dye}) * \epsilon_{prot} * D$$

**Without background and without dye correction:**

$$C = A_{280} * \epsilon_{prot} * D$$

C	Concentration in mg/ml
A <sub>280</sub>	Absorbance at 280 nm (10 mm path)
A <sub>320</sub>	Absorbance at 320 nm (10 mm path)
A <sub>max,dye</sub>	Absorbance value at the absorbance maximum of the dye (10 mm path)
ε <sub>prot</sub>	Extinction coefficient/protein factor in g*cm/l
cf <sub>dye</sub>	Dye-dependent correction factor at 280 nm
D	Dilution factor

- **Dye Concentration**

For dye-labeled proteins, the concentration of the dye is calculated using a modified form of the Beer-Lambert equation. For these calculations, the instrument considers the absorption maximum of the dye, and a dye-specific extinction coefficient (see Table 4). The dye concentration is calculated with or without background correction for are as follows:

**With background correction:**

$$C = \frac{((A_{max,dye} - A_{320}) * D)}{\epsilon_{dye} * 10^6}$$

**Without background correction:**

$$C = \frac{(A_{max,dye} * D)}{\epsilon_{dye} * 10^6}$$

C	Concentration
A <sub>max, dye</sub>	Absorbance at the max dye absorbance value (10 mm path)
A <sub>320</sub>	Absorbance at 320 nm (10 mm path)
ε <sub>dye</sub>	Extinction coefficient of dye in M <sup>-1</sup> *cm <sup>-1</sup>
D	Dilution factor

- **Degree of Labeling (DOL)**

DOL is the degree of labeling based on the average number of dye molecules coupled to a protein molecule. The degree of labeling can be determined from the absorption spectrum of the labeled antibody with or without background/dye correction as follows:

**With background and with dye correction:**

$$DOL = \frac{(A_{\max, \text{dye}} - A_{320}) * \epsilon_{\text{prot}}}{(A_{280} - A_{320} - cf_{\text{dye}} * (A_{\max, \text{dye}} - A_{320})) * \epsilon_{\text{dye}}}$$

**With background and w/o dye correction:**

$$DOL = \frac{(A_{\max, \text{dye}} - A_{320}) * \epsilon_{\text{prot}}}{(A_{280} - A_{320}) * \epsilon_{\text{dye}}}$$

**W/o background and with dye correction:**

$$DOL = \frac{A_{\max, \text{dye}} * \epsilon_{\text{prot}}}{(A_{280} - cf_{\text{dye}} * A_{\max, \text{dye}}) * \epsilon_{\text{dye}}}$$

**W/o background and w/o dye correction:**

$$DOL = \frac{A_{\max, \text{dye}} * \epsilon_{\text{prot}}}{A_{280} * \epsilon_{\text{dye}}}$$

DOL	Degree of labeling/dye per protein ratio
$A_{\max, \text{dye}}$	Absorbance value at the absorbance maximum of the dye (10 mm path)
$A_{280}$	Absorbance at 280 nm (10 mm path)
$A_{320}$	Absorbance at 320 nm (10 mm path)
$\epsilon_{\text{dye}}$	Extinction coefficient of dye in $M^{-1} * cm^{-1}$
$\epsilon_{\text{prot}}$	Extinction coefficient/protein factor in $g * cm/l$
$cf_{\text{dye}}$	Dye-dependent correction factor at 280 nm

**Table 4. Dye Types, Absorbance Max, Extinction coefficients, and dye-dependent correction factors**

NanoPhotometer Models	Dye Type	Absorbance maximum of Dye (nm)	Dye-dependent ext. coeff. $\epsilon_{\text{dye}}$ in $M^{-1} * cm^{-1}$	Dye-dependent correction factor at 280 nm $cf_{\text{dye}}$
NP80/N60/N50/C40	Alexa Fluor 350	346	19,000	0.19
NP80/N60/N50/C40	Alexa Fluor 405	401	34,500	0.70
NP80/N60/N50/C40	Alexa Fluor 488	494	71,000	0.11
NP80/N60/N50/C40	Alexa Fluor 532	530	81,000	0.09
NP80/N60/N50/C40	Alexa Fluor 546	554	104,000	0.12
NP80/N60/N50/C40	Alexa Fluor 555	555	150,000	0.08
NP80/N60/N50/C40	Alexa Fluor 568	577	91,300	0.46
NP80/N60/N50/C40	Alexa Fluor 594	590	73,000	0.56

NP80/N60/C40	Alexa Fluor 647	650	239,000	0.03
NP80/N60/C40	Alexa Fluor 680	679	184,000	0.05
NP80/N60/C40	Alexa Fluor 790	785	260,000	0.08
NP80/N60/N50/C40	Cy3	550	150,000	0.05
NP80/N60/C40	Cy5	649	250,000	0.05
NP80/N60/C40	DyLight 649	654	250,000	0.04
NP80/N60/N50/C40	DyLight 488	493	70,000	0.15
NP80/N60/N50/C40	FITC	495	68,000	0.30
NP80/N60/N50/C40	Pacific Blue	409	30,000	0.20
NP80/N60/N50/C40	Pacific Orange	397	24,500	0.60
NP80/N60/N50/C40	pHrodo Green	505	75,000	0.20
NP80/N60/N50/C40	pHrodo Red	560	65,000	0.12
NP80/N60/N50/C40	r-PE	566	200,000	0.18
NP80/N60/N50/C40	Texas Red	595	80,000	0.18

### ▪ Ratios

Protein samples e.g. from whole cell lysates may contain nucleic acids. To check the purity of the isolated protein, the 260/280 ratio is calculated to give an indication of the nucleic acid contamination. A pure protein preparation has an expected 260/280 ratios of ~ 0.57. If a ratio is detected out of the exseptable range an alert icon  is shown in the the results/table area. A click on the alert icon shows additional information.

The ratio is calculated with or without background correction according to if the background correction is activated during the measurements or not as follows:

#### Without background correction:

$$260/280 \text{ ratio} = \frac{A_{260}}{A_{280}}$$

#### With background correction:

$$260/280 \text{ ratio} = \frac{A_{260} - A_{320}}{A_{280} - A_{320}}$$

## PROTEIN ASSAYS

### METHOD OVERVIEW

Protein concentration may be measured using colorimetric assays, in which certain reagents are added to the protein solution to generate a colored product; either a protein-cupric ion chelate as in the Biuret, Lowry, BCA assays or a protein-dye complex as in the Bradford assay. In these colorimetric assays, the absorbance is measured in the visible range at the appropriate wavelength for each assay and compared against a standard curve prepared by serial dilution of a protein standard of known concentration. A linear regression analysis of the calibration standard data points is calculated by the NanoPhotometer. A correlation coefficient ( $R^2$ ) in the range of 0.95 to 1.00 indicates a good fit to a straight line.

- **Bradford Assay**

Method depends on quantifying the binding of a dye, Coomassie Brilliant Blue, to an unknown protein and comparing this binding to that of a standard curve prepared from a set of known protein of known concentrations at 595 nm. This standard is usually BSA (bovine serum albumin).

- **Biuret Assay**

Method depends on a reaction between cupric ions and peptide bonds in an alkali solution, resulting in the formation of a complex absorbing at 546 nm.

- **BCA Assay**

Method depends on a reaction between cupric ions and peptide bonds coupled with the detection of cuprous ions using bicinchoninic acid (BCA), giving an absorbance maximum at 562 nm. The BCA process is less sensitive to the presence of detergents used to solubilize membranes.

- **Lowry Assay**


Method is based on the Biuret reaction. Under alkaline conditions the divalent copper ion forms a complex with peptide bonds in which it is reduced to a monovalent ion. Monovalent copper ion and the radical groups of tyrosine, tryptophan, and cysteine react with Folin reagent to produce an unstable product that becomes reduced to molybdenum/tungsten blue. The bound reagent changes color from yellow to blue. This binding is compared with that obtained with a standard protein at 750 nm; this is usually BSA (bovine serum albumin).

**Note:** Detailed protocols are customarily supplied with these assay kits, and must be closely followed to ensure that accurate results are obtained.

## MEASUREMENT PROTOCOL

1. Select protein assays icon from home screen.  
To change between NanoVolume and cuvette application, use the Change to Cuvette/NanoVolume button below the parameter area (NP80/C40 only).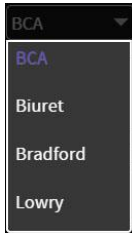

2. Select protein assay type from drop-down of the parameter area.  
Options are: BCA Assay, Bradford Assay, Lowry Assay, Biuret Assay

3. **For Cuvette Application:**  
Select the path length depending on the used cuvette.  
Options are: 1 mm, 2 mm, 5 mm and 10 mm  
**For NanoVolume Application:**  
Select the dilution depending on the sample concentration  
**Note:** There is no automatic path length setting available in this method.  
Select either a virtual dilution of 15 (path length 0.67 mm) or of 140 (path length 0.07 mm).


**For NanoVolume application with submicroliter cell (C40 only):**  
Select Lid depending on your sample concentration for concentration ranges of different lids (see also page 37 Submicroliter Cell Measurements Basics). Required sample volume for each lid is shown below the lid selection dropdown.

**Note:** A new Blank is recommended when changing the dilution lid.

4. If it is desired to heat the sample to 37°C use the toggle switch to turn on cell holder heater. When the cuvette holder has 37°C the color changes to green.

**Note:** Only available for cuvette applications (NP80 and C40).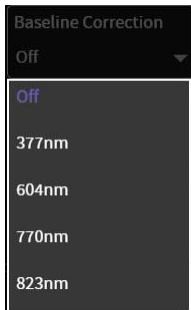


5. The baseline correction is disabled as the default. It is possible to choose between different wavelengths for the baseline correction.  
Options are: 377 nm, 604 nm, 770 nm (not N50) and 823 nm (not N50)

6. Select the curve fit type: Options are linear regression and zero regression (forces the straight line through the origin)
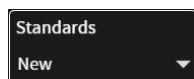

7. Standard new (no selection possible)

8. Select Unit

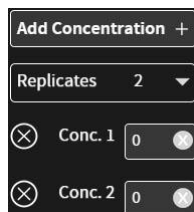9. **New:**  
Add up to 20 Concentrations and choose Replicates none, 2 or 3.

Measure a blank and depending on the replicate setting all concentrations. Results will be shown in the results area and if replicates are selected a mean value is shown. It is possible to exclude single measurements from the curve calculation by switching the toggle switch off.

**Note:** The standard curve cannot be altered any longer once the first sample is measured.

10. Once Standard curve is created or loaded it will be used for concentration calculations in the method. It might be necessary to do a blank measurement.

11. Apply sample and press the sample button to initiate the measurement.  
**Note:** Once the sample measurement is initiated it is no longer possible to do changes to the standard curve.

## CALCULATIONS

Protein concentration is determined using the standard curve by correlating absorbance values of samples with known concentration to calculate the concentration of the unknown sample. In order to maintain accuracy and precision please ensure that the  $R^2$  value of the standard curve is 0.95 or greater.

## KINETICS

## METHOD OVERVIEW

The Kinetics application is useful for: measuring initial rates of enzyme-catalyzed reactions, performing progress-curve analysis for complete reactions, calculating basic Michaelis-Menten parameters for single-substrate reactions, and measuring enzyme inhibition. Simple kinetics studies, where the change in absorbance is followed as a function of time at a fixed wavelength, can be readily performed with the NanoPhotometer. The rate of a chemical reaction can be


measured using spectrophotometric methods by studying the change in absorbance at a fixed wavelength as a function of time. These changes in absorbance reflect corresponding changes in the concentration of reactants or products as the reaction progresses. The rate of many chemical reactions can be markedly accelerated by the presence of catalysts, which remain chemically intact during the reaction. Catalysts in the case of biochemical reactions are generally represented by enzymes, which are specialized protein catalysts. However, a few examples of special reactions catalyzed by RNA molecules also exist. Studying the kinetics of a reaction can reveal important details of the catalytic mechanism involved in terms of sequence steps, transition state of reactants or nature of enzyme inhibitors.

## MEASUREMENT PROTOCOL

**Note:** If a kinetic is started from a control device via WiFi connection (tablet or smartphone) set the auto lock of the tablet or smartphone to never. Otherwise the kinetics will be interrupted when the smartphone or tablet is locked, because of losing the WiFi connection.

1. Select the Kinetics icon on the home screen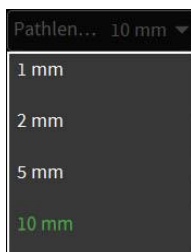

2. **For Cuvette Application:**

Select the path length depending on the used cuvette.

Options are: 1 mm, 2 mm, 5 mm and 10 mm

**For NanoVolume Application:**

Select the dilution depending on the sample concentration

**Note:** There is no automatic path length setting in this method. Select either a virtual dilution of 15 (path length 0.67 mm) or of 140 (path length 0.07 mm)

**For NanoVolume application with submicroliter cell (C40 only):**

Select Lid depending on your sample concentration for concentration ranges of different lids (see also page 37 Submicroliter Cell Measurements Basics). Required sample volume for each lid is shown below the lid selection dropdown.

**Note:** A new Blank is recommended when changing the dilution lid.


3. If it is desired to heat the sample to 37°C use the toggle switch to turn on cell holder heater. When the cuvette holder has 37°C the color changes to green.

**Note:** Only available for cuvette applications (NP80 and C40).

4. Default wavelength ( $\lambda$ ) is 340 nm but can be changed in the range of 200 – 900 nm, depending on the application.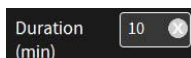


5. Enter the duration time in minutes over which measurements are taken. Possible range is 1 – 3000 minutes.

**Note:** A maximum of 500 samples is possible. Please consider by choosing the duration and interval time.

6. Enter the interval time between measurements in seconds. Possible interval times are 10 – 3,600 seconds, depending on the duration time.

**Note:** A maximum of 500 samples is possible. Please consider by choosing the duration and interval time.

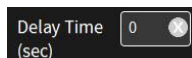7. Enter the delay time in seconds before the first measurement is taken. Possible delay time is between 0 – 3,600 seconds, depending on the duration time.

8. Insert cuvette with the reference sample and select blank button to initiate the measurement

9. Insert cuvette with the sample and select the sample button to initiate the measurements. Once the kinetic is started the Blank button turns to a Pause/Continue button and the Sample button to a Stop button.

**Note:** During the kinetics is running it is not possible to change the parameters, save data or delete data. Change of parameters is only possible before starting the kinetic. Save and delete data is only available after the kinetic is stopped.

## CALCULATIONS

All absorbance values are normalized to a 10 mm path.

$A_0$  = absorbance of start value (10 mm path)

$A_n$  = absorbance of actual value time n (10 mm path)

dA = absorbance of actual value – absorbance of start value

Slope = linear regression fit of all actual measurement points

Final A = absorbance of final value

$R^2 = R^2 = \sum_{i=1}^n \frac{(O_i - E_i)^2}{E_i}$  [ $O_i$  = observed slope value;  $E_i$  = expected slope value]

## OD600

### METHOD OVERVIEW

The growth of bacteria in liquid culture media is commonly monitored by measuring the optical density at 600 nm (OD600) in small samples taken from the cultures. OD600 measurements are typically used to determine the stage of growth of the bacterial culture, thereby ensuring that cells are harvested at an optimum point that corresponds to an appropriate density of live cells. Growth of bacterial cells typically progresses through a series of consecutive phases including: lag, log, stationary and decline (see Figure 1). In general, cells should be harvested towards the end of the log phase using the optical density of the samples to determine when this point has been reached. Since optical density in the case of OD600 measurements results from light scattering rather than light absorption, this value varies depending on the type of bacterial cells in the culture in terms of size and shape. Cells are routinely grown until the absorbance at 600 nm (known as OD 600) reaches approximately 0.4 prior to induction or harvesting. A linear

relationship exists between cell number (density) and OD 600 up to an absorbance value of 0.6, approximately.

As mentioned above, for turbid samples such as cell cultures, the absorbance measured is due to light scattering, and not the result of molecular absorption. Since the extent of scattering is affected by the optics of the system (distance between the cell holder and instrument exit slit, monochromator optics, slit geometry, etc.), different spectrophotometer types will tend to give different OD 600 readings for the same turbid sample. Therefore, if results from different spectrophotometers are to be compared, they must be normalized first using appropriate calibration curves.

A calibration curve can be constructed by comparing measured OD 600 to expected OD 600. Expected OD 600 is determined by counting cell number using an alternative technique (for example microscope slide method) and converting to OD 600 using the rule of thumb that  $1 \text{ OD } 600 = 5 \times 10^8 \text{ cells/ml}$  for *E. coli*.

The NanoPhotometer comes with a correction factor of 1 as default. To compare OD 600 values between different spectrophotometers, it is necessary to determine the constant deviation or ratio between the absorbance values for the same sample from each instrument and use this factor within the setting “correction factor” of your NanoPhotometer Software.

**Note:** The use of 10 mm path length disposable cuvettes is recommended for optical density measurements of cell culture solutions. The amount of cells is reflected in the reading and the likelihood of fluctuating amount of cells in a drop from sample to sample can be considered as extremely significant. It is therefore recommended to use cuvettes since the amount of error in a bigger volume is not as significant. The cuvette measurements provide a bigger average and therefore more reproducible readings. Also, to prevent the suspension settling too quickly and giving an OD reading that changes with time, glycerol should be added to the sample.


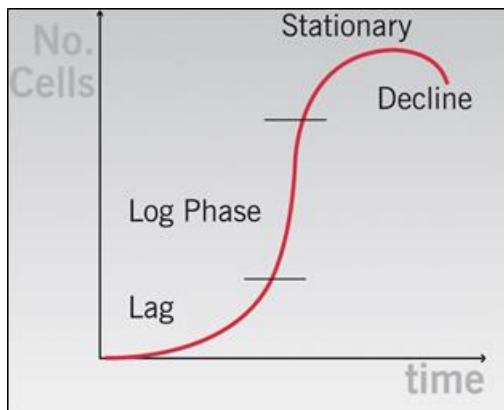
Figure 1 Bacterial growth curve

## MEASUREMENT PROTOCOL

1. Select the OD600 icon on the home screen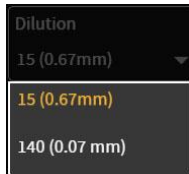

2. **For Cuvette Application:**

Select the path length depending on the used cuvette.  
Options are: 1 mm, 2 mm, 5 mm and 10 mm

**For NanoVolume Application:**

Select the dilution depending on the sample concentration

**Note:** There is no automatic path length setting in this method. Select either a virtual dilution of 15 (path length 0.67 mm) or of 140 (path length 0.07 mm)

**For NanoVolume application with submicroliter cell (C40 only):**

Select Lid depending on your sample concentration for concentration ranges of different lids (see also page 37 Submicroliter Cell Measurements Basics).

Required sample volume for each lid is shown below the lid selection dropdown.

**Note:** A new Blank is recommended when changing the dilution lid.

3. If it is desired to heat the sample to 37°C use the toggle switch to turn on cell holder heater. When the cuvette holder has 37°C the color changes to green.

**Note:** Only available for cuvette applications (NP80 and C40).


4. Default wavelength is 600 nm but can be changed in the range of 200 – 900 nm, depending on the application.

5. Toggle switch cells/ml is default disabled. Enable cells/ml to get the cells/ml calculated. Enter the cell specific factor and multiplier (e.g. 1 OD600 = 5 x 10<sup>8</sup> cells/ml)

6. Enter the correction factor to compensate for different optical configurations between this and other instruments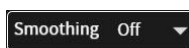

7. Option to smooth the graph with different boxcars. off, 1 = boxcar 11 (default), 2 = boxcar 21 and 3 = boxcar 61

8. Option to set/calculate a dilution factor for manual diluted samples.

9. Insert cuvette with the reference sample and select blank button to initiate the measurement

10. Insert cuvette with the sample and select the sample button to initiate the measurement

## CALCULATIONS

$$OD600 = A_{600} * D * cf$$

OD600	Optical density at 600 nm
$A_{600}$	Absorbance at 600 nm (10 mm path)
D	Dilution factor
cf	Correction factor for spectrophotometer

$$\text{Cells/ml} = A_{600} * D * cf * \text{multiplier}$$

$A_{600}$	Absorbance at 600 nm (10 mm path)
D	Dilution factor
cf	Correction factor for spectrophotometer
multiplier	Multiplier of sample

## MORE APPS

The More Apps icon located on the home screen opens another menu screen with access to icons for additional applications available on the NanoPhotometer. The applications featured in this menu include: wavelength, concentration, wavescan, absorbance/ratio, standard curve and custom applications.

## MORE APPS: WAVELENGTH

### METHOD OVERVIEW

In the wavelength application it is possible to measure simple absorbance (A) and % transmittance (%Trans. / only in cuvette mode) of a sample at specific wavelengths. It is possible to add up to 20 different wavelengths

### MEASUREMENT PROTOCOL

1. Select the More Apps icon  from the Home screen and then the Wavelength icon from the More Apps screen


2. **For NanoVolume application:**  
Select the dilution depending on the sample concentration  
**Note:** There is no automatic path length setting in this method. Select either a virtual dilution of 15 (path length 0.67 mm) or of 140 (path length 0.07 mm)  
**For cuvette application:**  
Select the path length depending on the used cuvette.

Options are: 1 mm, 2 mm, 5 mm and 10 mm

**For NanoVolume application with submicroliter cell (C40 only):**


Select Lid depending on your sample concentration for concentration ranges of different lids (see also page 37 Submicroliter Cell Measurements Basics). Required sample volume for each lid is shown below the lid selection dropdown.

**Note:** A new Blank is recommended when changing the dilution lid.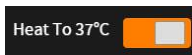


- If it is desired to heat the sample to 37°C use the toggle switch to turn on cell holder heater. When the cuvette holder has 37°C the color changes to green.


**Note:** Only available for cuvette applications (NP80 and C40).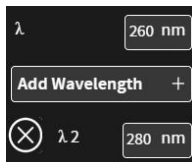


- Enter desired wavelength ( $\lambda$ ) to be measured. It is possible to measure up to 20 wavelengths simultaneously. More wavelength ( $\lambda$ ) options are added by selecting the wavelength button. An added wavelength can be deleted with  $\otimes$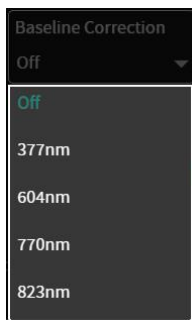

- Baseline correction is default off and can be enabled by selecting the different values for the correction in the drop-down menu: 377 nm, 604 nm, 770 nm (N/A N50) and 823 nm (N/A N50)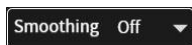

- Option to smooth the graph with different boxcars. off, 1 = boxcar 11 (default), 2 = boxcar 21 and 3 = boxcar 61

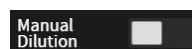- Option to set/calculate a dilution factor for manual diluted samples.

- Apply the blank ddH<sub>2</sub>O or buffer to the illuminated sample window for the reference measurement and select blank to initiate the reading.

**Note:** The illumination of the sample window can be switched off in the preferences.

- Use a lint-free laboratory wipe to clean both the sample window and mirror prior to applying the next sample.

**Note:** It could be helpful to apply the blank a second time and read it as a sample to ensure a proper blank.

- Apply sample to the sample window and press the sample button to initiate

the measurement.

## CALCULATIONS

### ▪ Absorbance Calculation:

Absorbance is formally defined as the decimal logarithm (base 10) of the reciprocal of transmittance:

$$A = \log\left(\frac{1}{T}\right) = -\log T$$

$$T = 10^{(-A)}$$

**Note:** Corresponding absorbance value e.g. Absorbance value ( $\lambda = 230$ ) etc. normalized to 10 mm path length

### ▪ %Transmittance Calculations (cuvette mode only)

In the wavelength application it is possible to measure the absorbance (A) and % transmittance (%T) on a sample with respect to a reference at a specific wavelength. Transmittance is the ratio of light intensity remaining after it has passed through the sample ( $I$ ) to the initial incident light intensity ( $I_0$ ):

$$T = \frac{I}{I_0}$$

$$\%T = \frac{I}{I_0} \times 100$$

## MORE APPS: WAVESCAN

### METHOD OVERVIEW

Using the wavescan application it is possible to obtain the complete spectral scan for a defined wavelength range between 200 - 900 nm (NP80/N60/C40) or from 200 - 650 nm (N50).

### MEASUREMENT PROTOCOL

1. Select the More Apps icon  from the Home screen and then the Wavescan icon from the More Apps screen

2. **For NanoVolume application:**

Select the dilution depending on the sample concentration

**Note:** There is no automatic path length setting in this method. Select either a virtual dilution of 15 (path length 0.67 mm) or of 140 (path length 0.07 mm)

**For cuvette application:**

Select the path length depending on the used cuvette.

Options are: 1 mm, 2 mm, 5 mm and 10 mm

**For NanoVolume application with submicroliter cell (C40 only):**

Select Lid depending on your sample concentration for concentration ranges of different lids (see also page 37 Submicroliter Cell Measurements Basics).

Required sample volume for each lid is shown below the lid selection dropdown.

**Note:** A new Blank is recommended when changing the dilution lid.


3. If it is desired to heat the sample to 37°C use the toggle switch to turn on cell holder heater. When the cuvette holder has 37°C the color changes to green.

**Note:** Only available for cuvette applications (NP80 and C40).


4. Set desired wavelength range (Start and End Wavelength) to be scanned for the samples.

**Note:** If samples with a different wavelength range are selected, the graphs are shown on the full scan range of 200 – 900 nm (N50: 200 - 650 nm)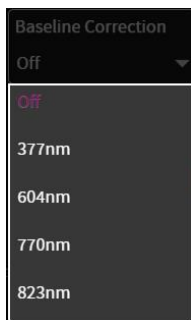

5. Baseline correction is off as the default and can be enabled by selecting the different values for the correction in the drop-down menu: 377 nm, 604 nm, 770 nm (N/A N50) and 823 nm (N/A N50)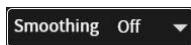

6. Option to smooth the graph with different boxcars. off, 1 = boxcar 11 (default), 2 = boxcar 21 and 3 = boxcar 61


7. Option to set/calculate a dilution factor for manual diluted samples.

8. Apply the blank ddH<sub>2</sub>O or buffer to the illuminated sample window for the reference measurement and select blank to initiate the reading.

**Note:** The illumination of the sample window can be switched off in the preferences.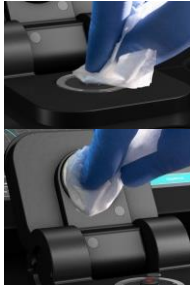

- Use a lint-free laboratory wipe to clean both the sample window and mirror prior to applying the next sample.  
**Note:** It could be helpful to apply the blank a second time and read it as a sample to ensure a proper blank.

- Apply sample to the sample window and press the sample button to initiate the measurement.

## CALCULATIONS

No calculations necessary, values are reported based on 10 mm path length.

In the results are the prominent peaks shown with wavelength and absorbance value. For cuvette measurements it is possible to change to the %Transmittance mode. If a peak of interest is not shown in the results the peak can be added by clicking on the graph and added to the results by tapping on the Add Peak button in the pop-up window:

Add peak	
Sample Name	Sample22
Wavelength	259
Absorbance	0.771

## MORE APPS: ABSORBANCE RATIO

### METHOD OVERVIEW

In this mode, it is possible to determine simple absorbance ratios for a given sample by measuring the absorbance at two wavelengths specified in the parameters of the method relative to a blank.

### MEASUREMENT PROTOCOL

- Select the More Apps icon  from the Home screen and then the Absorbance/Ratio icon from the More Apps screen

2. **For NanoVolume application:**

Select the dilution depending on the sample concentration

**Note:** There is no automatic path length setting in this method. Select either a virtual dilution of 15 (path length 0.67 mm) or of 140 (path length 0.07 mm)

**For cuvette application:**

Select the path length depending on the used cuvette.

Options are: 1 mm, 2 mm, 5 mm and 10 mm

**For NanoVolume application with submicroliter cell (C40 only):**


Select Lid depending on your sample concentration for concentration ranges of different lids (see also page 37 Submicroliter Cell Measurements Basics). Required sample volume for each lid is shown below the lid selection dropdown.

**Note:** A new Blank is recommended when changing the dilution lid.


3. If it is desired to heat the sample to 37°C use the toggle switch to turn on cell holder heater. When the cuvette holder has 37°C the color changes to green.

**Note:** Only available for cuvette applications (NP80 and C40).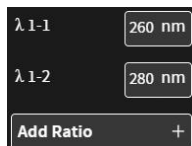

4. Enter desired wavelengths ( $\lambda$  1-1 and  $\lambda$  1-2) for ratio calculation. It is possible to measure up to 20 absorbance/ratios simultaneously. More wavelengths for ratio calculation can be added by selecting the Add Ratio button. The added ratios can be deleted with  $\otimes$

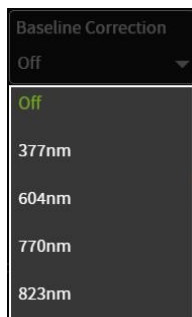5. Baseline correction is off as the default and can be enabled by selecting the different values for the correction in the drop-down menu: 377 nm, 604 nm, 770 nm (N/A N50) and 823 nm (N/A N50)

6. Option to smooth the graph with different boxcars. off, 1 = boxcar 11 (default), 2 = boxcar 21 and 3 = boxcar 61

7. Option to set/calculate a dilution factor for manual diluted samples.


8. Apply the blank ddH2O or buffer to the illuminated sample window for the reference measurement and select blank to initiate the reading.

**Note:** The illumination of the sample window can be switched off in the preferences.

9. Use a lint-free laboratory wipe to clean both the sample window and mirror prior to applying the next sample.

**Note:** It could be helpful to apply the blank a second time and read it as a

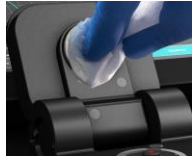sample to ensure a proper blank.

10. Apply sample to the sample window and press the sample button to initiate the measurement.

## CALCULATIONS

The absorbance ratio is calculated from the two path lengths specified by the user in the parameters.

$$\lambda_1 : \lambda_2 = \frac{\lambda_1}{\lambda_2}$$

$\lambda_1 : \lambda_2 =$  Absorbance Ratio

$\lambda_1 =$  Absorbance 1 corresponding absorbance value 1 selected normalized to 10 mm path

$\lambda_2 =$  Absorbance 2 corresponding absorbance value 2 selected normalized to 10 mm path

## MORE APPS: CONCENTRATION

### METHOD OVERVIEW

In this mode, concentration can be calculated for a sample by determining the absorbance at a specific wavelength relative to a reference. The concentration is then obtained by multiplying the measured absorbance by a specific factor. This factor may be known in advance and entered by the user, or it may be calculated by the instrument by measuring a set of standard (standard curve method) with known concentrations to create a standard curve.

### MEASUREMENTS PROTOCOL

1. Select the More Apps icon  from the Home screen and then the Concentration icon from the More Apps screen.

2. **For NanoVolume application:**  
 Select the dilution depending on the sample concentration  
**Note:** There is no automatic path length setting in this method. Select either a virtual dilution of 15 (path length 0.67 mm) or of 140 (path length 0.07 mm)  
**For cuvette application:**  
 Select the path length depending on the used cuvette.  
 Options are: 1 mm, 2 mm, 5 mm and 10 mm  
**For NanoVolume application with submicroliter cell (C40 only):**

Select Lid depending on your sample concentration for concentration ranges of different lids (see also page 37 Submicroliter Cell Measurements Basics). Required sample volume for each lid is shown below the lid selection dropdown.

**Note:** A new Blank is recommended when changing the dilution lid.

- If it is desired to heat the sample to 37°C use the toggle switch to turn on cell holder heater. When the cuvette holder has 37°C the color changes to green.

**Note:** Only available for cuvette applications (NP80 and C40).

- Default wavelength is 260 nm but can be changed, depending on the application.

- Enter a factor for concentration calculation

- Baseline correction is off as the default and can be enabled by selecting the different values for the correction in the drop-down menu: 377 nm, 604 nm, 770 nm (N/A N50) and 823 nm (N/A N50)

- Option to smooth the graph with different boxcars. off, 1 = boxcar 11 (default), 2 = boxcar 21 and 3 = boxcar 61


- Option to set/calculate a dilution factor for manual diluted samples.

- Apply the blank ddH2O or buffer to the illuminated sample window for the reference measurement and select blank to initiate the reading.

**Note:** The illumination of the sample window can be switched off in the preferences.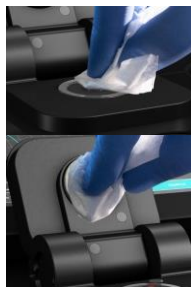

- Use a lint-free laboratory wipe to clean both the sample window and mirror prior to applying the next sample.

**Note:** It could be helpful to apply the blank a second time and read it as a sample to ensure a proper blank.

- Apply sample to the sample window and press the sample button to initiate the measurement.

## CALCULATIONS

In this method, the concentration of the sample is calculated based on the Beer-Lambert law given the user specified wavelength of interest and user defined extinction coefficient. The equations for calculating concentration without background correction are as follows:

**Without background correction:**

$$C = A_n * \epsilon * D$$

C	Concentration (ng/μl)
$A_n$	Absorbance at user specified path length n (10 mm path)
D	Dilution factor
$\epsilon$	extinction coefficient/factor

## MORE APPS: STANDARD CURVE


### METHOD OVERVIEW

The construction of a calibration curve from multiple standards of known concentrations can be created and stored on the NanoPhotometer. The standard curve can be used to quantify samples of the same type with unknown concentrations. This application provides an extremely useful tool with which to integrate, expedite and simplify the measurement and calculations involved in determining the concentration of analytes in unknown samples. If a zero concentration standard is required, include it in the number of standards to be entered using a reagent blank and entering 0.00 for concentration.

### MEASUREMENTS PROTOCOL

1. Select the More Apps icon  from the Home screen and then the Standard Curve icon from the More Apps screen.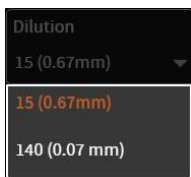

2. **For NanoVolume application:**  
Select the dilution depending on the sample concentration  
**Note:** There is no automatic path length setting in this method. Select either a virtual dilution of 15 (path length 0.67 mm) or of 140 (path length 0.07 mm)  
**For cuvette application:**  
Select the path length depending on the used cuvette.  
**For NanoVolume application with submicroliter cell (C40 only):**  
Select Lid depending on your sample concentration for concentration ranges of different lids (see also page 37 Submicroliter Cell Measurements Basics). Required sample volume for each lid is shown below the lid selection dropdown.

**Note:** A new Blank is recommended when changing the dilution lid.

3. If it is desired to heat the sample to 37°C use the toggle switch to turn on cell holder heater. When the cuvette holder has 37°C the color changes to green.


**Note:** Only available for cuvette applications (NP80 and C40).


4. The baseline correction is disabled as the default. It is possible to choose between different wavelengths for the baseline correction. Options are: 377 nm, 604 nm, 770 nm (N/A N50) and 823 nm (N/A N50)


5. Select the curve fit type: Options are linear regression and zero regression (forces the straight line through the origin)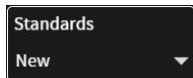

6. Standard new no selection possible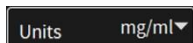

12. Select Unit

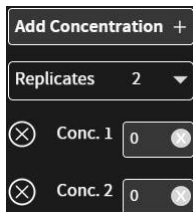**New:**

Add up to 20 Concentrations and choose Replicates none, 2 or 3.

Measure a blank and depending on the replicate setting all concentrations. Absorbances will be shown in the results area and if replicates are selected a mean value is shown. It is possible to exclude single measurements from the curve calculation by switching the toggle switch off.

**Note:** The standard curve cannot be altered any longer once the first sample is measured.

7. Once Standard curve is created or loaded it will be used for concentration calculations in the method. It might be necessary to do a blank measurement.

8. Apply sample and press the sample button to initiate the measurement. **Note:** Once the sample measurement is initiated it is not possible to do changes to the standard curve.

## CALCULATIONS

Concentration is determined via the absorbance values provided by the standard curve based on the curve fit selection including the following options: linear regression or zero regression.


## CUSTOM APPS

There is an option for designing customer specific Custom Apps which can be loaded to the NanoPhotometer. For more information about designing custom applications to suit individual research needs please contact Implen directly for assistance.

## STORED RESULTS

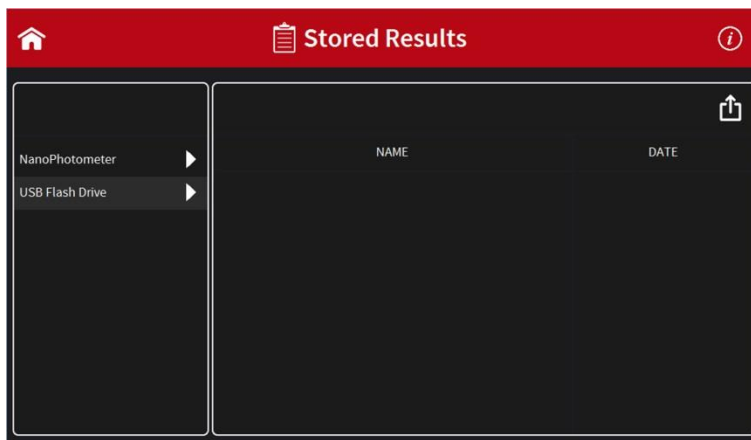The Stored Results icon opens the directories of folders containing files of results that have been previously saved.

On the left side of the screen all available directories/storages are shown: NanoPhotometer, Control Device and/or USB flash drive. A click on a storage shows on the left side the subfolders of the storage and on the right side the files. On the right side of the screen all saved result files of the selected folder are shown and can be opened by a long or double click. Folders can be deleted, renamed, moved or copied to by clicking on the  icon. It is also possible to delete, rename, move or copy files by clicking on the  icon.

**Note:** PDF and Excel files cannot be opened on the NanoPhotometer. The files need to be transferred to a computer or device where Excel or a PDF reader is installed.

**Note:** Control device is only available on coputer, tablets and smartphones not on the NanoPhotometer version of the software.


For data transfer via Ethernet or WiFi see page 40 Data Transer.

## STORED METHODS

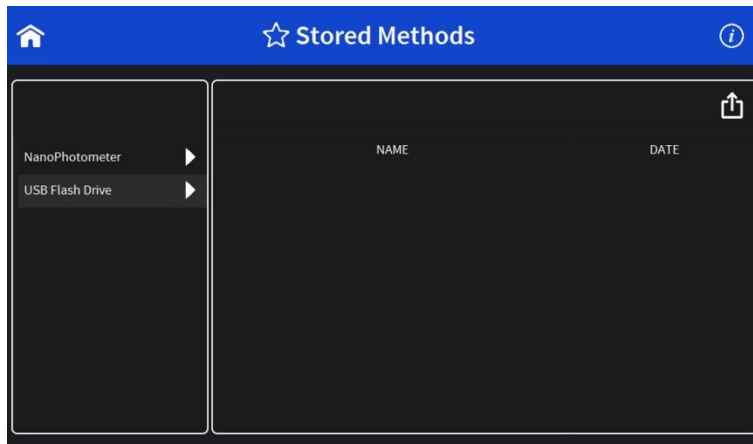The stored methods icon opens the directories of folders containing methods stored by the user.

On the left side of the screen all available directories/storages are shown: NanoPhotometer and/or USB flash drive. A click on a storage shows on the left side the subfolders of the storage and on the right side the method files. On the right side of the screen all saved method files of the selected folder are shown and can be opened by a long or double click.

New folders can be created by clicking on . Folders can be deleted, renamed, moved or copied to by clicking on the  icon. It is also possible to delete, rename, move or copy methods by clicking on the  icon.

## 5. PREFERENCES

System preferences can be set by selecting preferences () on the home screen. The preferences menu includes: general, dyes, warning messages and network. The menu item which is selected will have its preference options listed in the window to the right.

**Note:** Preferences are not available for smartphone < 7" screen versions.

### GENERAL

Selecting General in the Preferences menu opens a window to the right of the preferences menu with the following options: Display (built-in display only), About, WiFi Access Point and Illumination Sample Window (NP80/N60/N50) or Enable NanoVolume (C40).

### DATE AND TIME

Within the Date and Time it is possible to set the actual date and time. Change the date and/or time by clicking on the date/time field. The selected time is shown below the date and time field.

To change the time it is necessary to reboot the NanoPhotometer by clicking on the Set and Reboot button.

## DISPLAY

The brightness of the built-in screen can be adjusted under display.

**Note:** Display preferences are only available for instruments with integrated touchscreen.

## ABOUT

In About the following information of the NanoPhotometer are shown: NanoPhotometer Version, Serial Number, IP Address, Hardware Version, Firmware Version, Software Version, Time & Date of Initialization Test and Status of Selftest.

## WiFi ACCESS POINT

Toggle Switch to en-/disable the WiFi access point.

## ILLUMINATION SAMPLE WINDOW

Toggle switch to switch on/off the illumination of the sample window (for NP80/N60/N50 only)


**Note:** Only available for NP80/N60/N50 models.

## ENABLE NANOVOLUME

Toggle switch to enable the submicroliter cell option in the NanoPhotometer C40.

## DYES


There is a list of available dye-labels for both nucleic acid dye and protein dye methods. To toggle between the list click on the Nucleic Acid/Protein buttons in the header.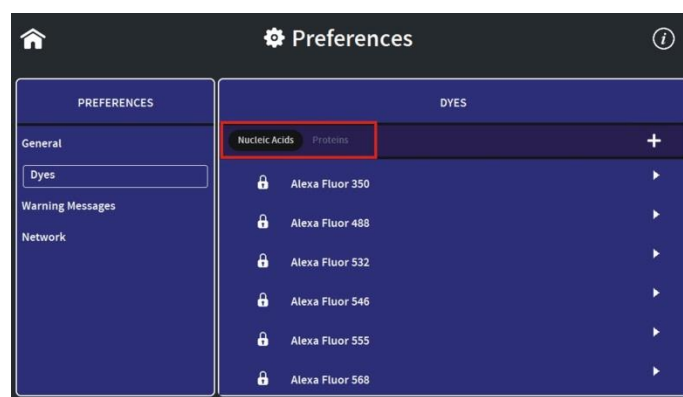

Each dye has either an (🔒) icon in front of the dye name indicating that the dye is locked and cannot be changed or deleted. If the dye has a delete symbol (🗑️), there is an option to delete the dye which is not locked or pre-programmed in the NanoPhotometer software.


Selecting a dye name opens a new screen with the dye information: dye name, absorbance maximum dye (nm), dye-dependent extinction coefficient  $\epsilon_{\text{dye}}$  ( $\text{M}^{-1} * \text{cm}^{-1}$ ), and dye-dependent

correction factor as well as the option to show the dye in the drop-down of the application (Nucleic Acid or Protein UV).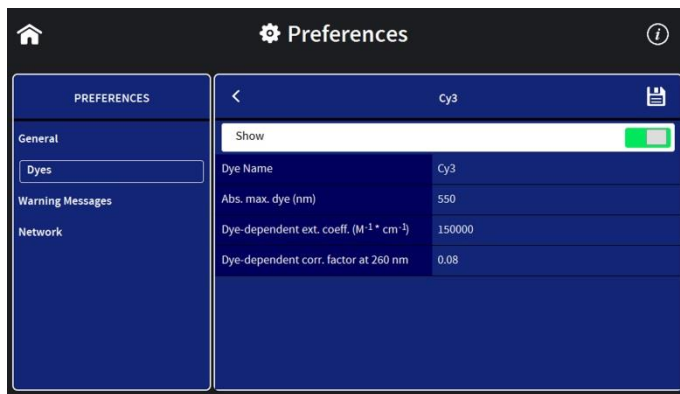

**Note:** It is not possible to delete a dye of the factory list; custom dyes can be deleted if they are not locked.

It is possible to add a new dye to the list by selecting the **+** button to add a new dye. A window will open where it is possible to enter the: dye name, dye absorbance maximum (nm), dye-dependent extinction coefficient  $\epsilon_{dye}$  ( $M^{-1} \cdot cm^{-1}$ ), and dye-dependent correction factor. There is a toggle switch available to lock the dye to prevent deleting a dye from the dye list accidentally.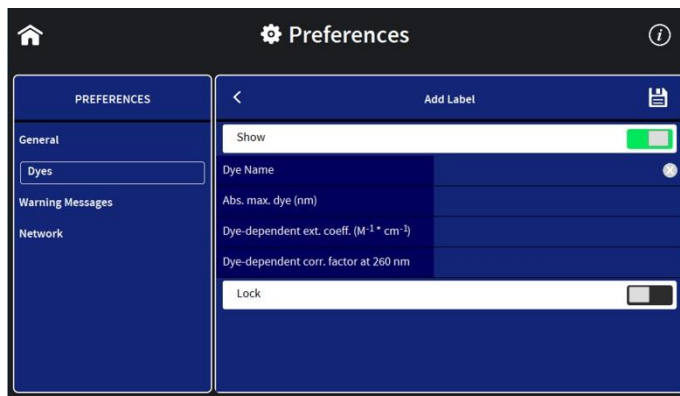


## WARNING MESSAGES

### BLANK CONTROL


Toggle switch to switch on/off the Blank Control™ of the NanoPhotometer.

**Note:** The Blank Control™ is available for all NanoVolume methods (for NP80/N60/N50).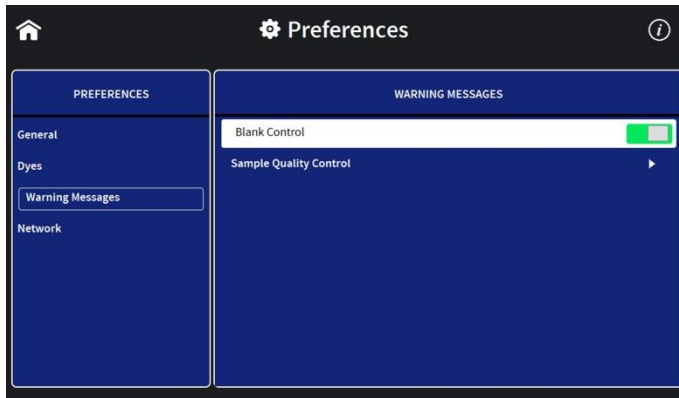


## SAMPLE QUALITY CONTROL

It is possible to change the upper and the lower limit of the ratio alert warning messages. Default values for Nucleic Acids are for 260/230 ratio 1.8 A - 3 A and for 260/280 ratio 1.65 A - 2.5 A. The upper limit for the Protein UV 260/280 ratio is 0.7 A.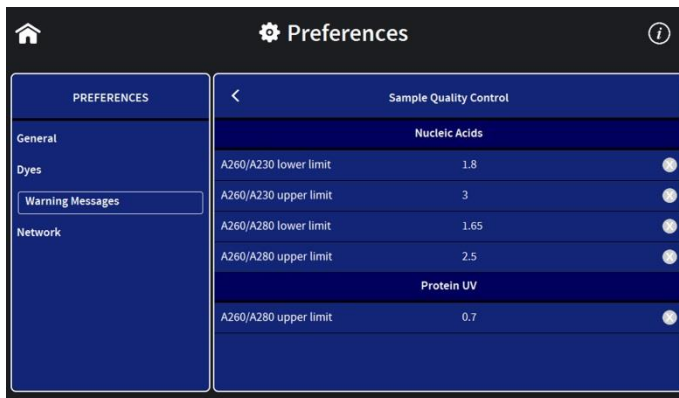

## NETWORK

In Network preferences it is possible to choose between a Dynamic Host Configuration Protocol (DHCP) and a static network configuration. As default the DHCP protocol is selected. Connect the NanoPhotometer with the Ethernet and the IP is automatically set during the start up of the NanoPhotometer. If none IP is shown click on the save button.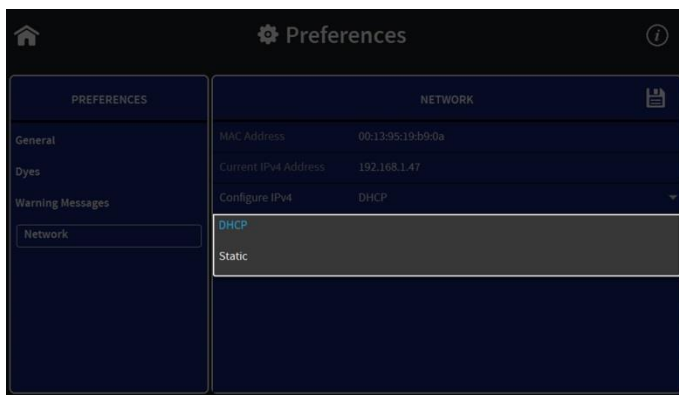

For static IP configuration select Static in the Configure IPv4 dropdown and enter the IPv4 address, netmask and gateway and confirm with the save button.

## 6. TROUBLESHOOTING

### SELF-CALIBRATION TEST

The NanoPhotometer self-calibration test is performed automatically every time the instrument is powered on. If the instrument passes the self-calibration test the home screen is shown. If the instrument does not pass the test, a window with a message will appear along with the message explaining the reason for the failed test along with the recommended solution. If the ok button is selected, the window is closed and the home screen will be shown with an error message at the top. The error message is also shown on top of all other screen and all performed measurements including the saved files of those measurements. If the self-calibration test fails please contact the Implen Support Team.


## 7. ASSISTANCE

The Assistance menu includes: support, report a problem (only available for tablet and computer versions), software maintenance and legal as functions to help with any technical issues or questions that may arise with the NanoPhotometer.

**Note:** Assistance function is not available for smartphone < 7" screen versions.

### SUPPORT

Selecting support on the left side assistance menu will show the available options for contacting Implen.

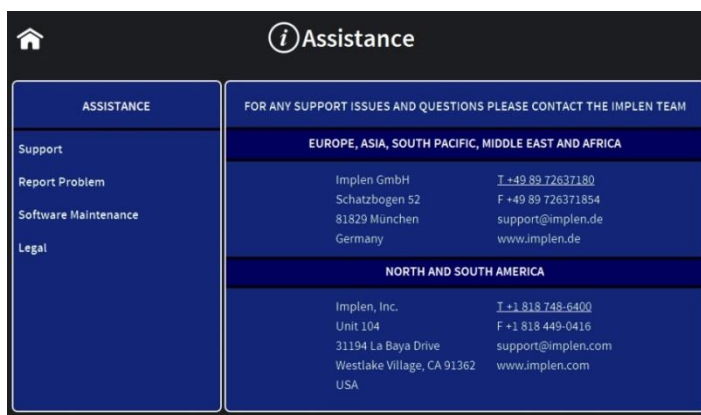### REPORT PROBLEM

The function to report a problem is available on computer and tablet versions only. When report a problem is highlighted in the left side assistance menu, the right side will have a form to fill out

with information including: first name, last name, phone number, email, and country. A drop-down menu provides the option to select the problem type and includes the following selections: error message, software, firmware, measurements, and other. It is also possible to enter a question or comment at the end of the form. Once the form is completed and the send button is selected a message will be sent directly to Implen and the appropriate support person will contact as soon as possible to provide further support.

**Note:** The report a problem function is only available for computer and tablet versions.

## SOFTWARE MAINTENANCE

### RESET

There is an option to reset the instrument to factory settings. By selecting the reset button a window will open that says “Reset the NanoPhotometer?” Selecting the cancel button will close the window without changing the settings and selecting reset will open a window that will ask again “Reset the NanoPhotometer to factory settings? All data, stored methods and settings will be lost.” If it is confirmed the factory settings will be restored.

**Note:** All stored methods, settings and data on the NanoPhotometer will be deleted if the reset option is executed.

### SOFTWARE UPDATE

Download the software update file from the Implen homepage: [www.implen.de/downloads/](http://www.implen.de/downloads/) and copy it to a USB flash drive.

**Note:** Make sure that the sample name is NPOS.bin without any additions or extensions like from multiple downloads, otherwise rename file to be NPOS.bin exactly.

**Note:** Save all data on the NanoPhotometer before updating.

**Note:** Make sure that the NanoPhotometer is connected to power and that the power connection is not interrupted during the update.

It is recommended to do the update via the built-in touchscreen. If the touchscreen is not available it is also possible to do the update via computer or tablet. In this case it is necessary to relaunch the NanoPhotometer software after the reboot of the NanoPhotometer. For tablets the WiFi connection needs to be reconnected.

**Note:** Update always first the firmware of the NanoPhotometer and then the client software (PC/MAC, smartphone, tablet).

#### Update procedure:

1. Copy the NPOS.bin file to a USB flash drive into the root folder
2. Insert in the USB flash drive to the USB port of the NanoPhotometer
3. Select Assistance/Software Maintenance
4. Click on “Update” and wait until the NanoPhotometer reboots

**Note:** Please update also in a second step the client software (PC/MAC, tablet, smartphone).

## CREATE LOG FILE

To create a log file insert a USB flash drive to the USB port of the NanoPhotometer. Click on “Create Log File”. The log (NPOS.log) file will be saved on the root folder of the USB flash drive.

## INSTALL PRINTER DRIVER

Printer driver installation is only necessary if the NanoPhotometer runs with a software version lower than 1.1.11147.

1. Switch NanoPhotometer on / home screen
2. Update to newest firmware version

### **Update Process:**

- a. Copy NPOS.bin to a USB flash drive into the root folder
- b. Insert the USB flash drive into your NanoPhotometer
- c. Select: “Assistance / Software Maintenance”
- d. Click on “Update”
- e. Wait until the NanoPhotometer reboots

**Note:** Update always first the firmware on the NanoPhotometer and in a second step the client software (PC/MAC, tablet, smartphone).

3. Go to “Assistance / Software Maintenance”
4. Start driver installation by a click to the “Install Driver for Printers” button
5. Wait until the following message is shown: “Installed driver for printer successfully.”
6. Go to home screen
7. Connect Dymo/HP printer via USB cable
8. Wait for 30 seconds
9. Dymo/HP printer is ready to use

**Note:** Connect the printer after starting the NanoPhotometer or in the home screen. It is not possible to operate more than one printer at a time.

## LEAGAL

The NPOS software is Copyright of Implen Inc. and its affiliates. The NPOS software includes some open source software components under the licenses listed in “Legal”.

## CONTACT IMPLEN

There is an option to contact Implen for instruments connected to the internet. For any support issues and questions please contact the Implen team directly:

**Europe, Asia, South Pacific, Middle East and Africa**

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## 8. MAINTENANCE

### MAINTENANCE FREE TECHNOLOGY

The NanoPhotometer technology is maintenance-free. Regular maintenance and calibration is not necessary.

For facilities that are working according to national as well as international guidelines and standards including: Good Laboratory Practice (GLP), Good Manufacturing Practice (GMP), or ISO9000-9004; the proper performance of the spectrophotometer has to be tested and proven on a regular basis with individually set intervals. Implen provides certified NanoPhotometer secondary standards as an optional accessory. These NanoPhotometer Didymium glass filters and standard solution sets are suitable for the control and documentation of the wavelength accuracy and the photometric accuracy of your system. IQ/OQ documentation is also available. Please contact your local Implen office or an authorized Implen partner for further information.

Support agreements that help to fulfill the demands of regulatory guidelines concerning GLP/GMP include: calibration certification using filters traceable to international standards (during production and quality control), certified engineers and calibrated test equipment,

approved to ISO 9001 standard, automatic self-diagnostic calibration test during start of the NanoPhotometer, result is documented in each data output file, and possibility to save a Implen Document Source (IDS) file (no data manipulation possible).

## REPLACEMENT PARTS

### ▪ Lamp Replacement

The NanoPhotometer is equipped with a xenon flash lamp and should not need replacement until several years of use. In the unlikely event the lamp does need to be replaced, this should be done by the manufacturer or a certified service engineer from your supplier.

### ▪ Battery Replacement

The optional battery pack can only be assembled or replaced by the manufacturer or a certified service engineer from your supplier.

**Note:** Danger of explosion if battery is incorrectly replaced. Replacement only with the same or equivalent type recommended by the manufacturer: Replacement needs to be done by a certified service engineer

### ▪ Touchscreen Replacement

The optional touchscreen can only be assembled or replaced by the manufacturer or a certified service engineer from your supplier.

## CLEANING AND GENERAL CARE

Switch off the NanoPhotometer and disconnect the power cord prior to external cleaning. Use a soft wet cloth or dry microfiber cloth to clean all external surfaces. A mild liquid detergent may be used to remove stubborn marks.

Approved disinfectant solutions include: Apesin disinfection spray (Tana Chemi GmbH), Incidin Liquid & Inciddin Foam (Ecolab), and Lysoformin Spezial (Lysoform Dr. Hans Roseman GmbH).

**Note:** Observe all necessary precautions if dealing with hazardous samples or solvents.

## MESSAGES

The software gives three different kind of messages confirmation (green), warning (yellow) and alert (red) messages.

Confirmation messages are shown e.g. after the initialization test has passed successfully or if files or folders are copied or moved successfully. These messages are self-explaining and not listed in the user manual.

**IMPORTANT WARNING MESSAGES:**

Message Text	Explanation/Solution
Air bubble, lint residue or bad sample. Please reapply sample.	Air bubble recognition is on and has detected either an air bubble, lint residue or a bad sample like e.g. a turbid sample. Check sample, clean the pedestal and mirror in the lid thoroughly and reapply sample carefully. To avoid air bubbles apply sample by reverse pipetting.
Another control device is currently connected.	Message is shown on a control device when another control device is currently connected with an open measurement session. Please close method on control device.
Battery is low. Please connect to power!	The message pops up twice. Between the first and the second message is at least one hour left to use the NanoPhotometer under standard conditions. After the second warning the instrument is automatically shutting down within 10 to 15 seconds.
Concentration too high. Change to lid10/lid50/lid100/lid250	The used sample is too high concentrated for submicroliter cell measurements with the used lid. Change to the recommended lid. A new blank is recommended when changing the dilution lid.
Concentration too high. Dilute sample.	The used sample is too high concentrated for submicroliter cell measurements Lid 250 with the C40. Sample is above the detection limit of the instrument. Dilute sample manually to measure within the detection range of the instrument.
Concentration too low.	The used sample is too low concentrated for submicroliter cell measurements Lid 5 with the C40. Sample is below the detection limit of the instrument. Only cuvette measurement may be possible
Concentration too low. Change to lid5/lid10/lid50/lid100	The used sample is too low concentrated for submicroliter cell measurements with the used lid. Change to the recommended lid. A new blank is recommended when changing the dilution lid.
High absorbance at 250-280nm, 280-340nm, 340-400nm, 400-475nm, 475-550nm, 550-625nm, 625-700nm. Bad Blank or insufficient cleaning.	The warning message is shown for when during a blank measurement (NanoVolume) an absorbance of $\geq 2.3$ (10 mm path) is detected. The warning message shows the range(s) at what range the absorbance is appearing. A high absorbance in a blank measurement can have two reasons: either the blank solution/buffer has an absorbance in this wavelength range or the pedestal was not cleaned properly after the last reading. Clean the pedestal and the mirror in the lid thoroughly and check the blank solution for absorbance (Use water for blank and measure the blank buffer as sample).
Implen Folder is Empty	Is shown on control devices like tablet and smartphone when no data are saved in the control device folder. For Android tablet and smartphone this folder is always empty. Files are saved in the download folder of the Android tablet smartphone directory.

Instrument connected to control device! Do you really want to intercept?	Message is shown on the NanoPhotometer and a method is open on another control device. Click on OK and the measurements session is interrupted and the data are saved in the auto save folder.
Lint residue or bad sample. Please reapply sample.	Air bubble recognition is off. Software has detected either a lint residue or a bad sample like e.g. a turbid sample. Check sample, clean the pedestal and mirror in the lid thoroughly and reapply sample carefully.
Maximum absorbance level at specified wavelength reached. Calculations may lead to low/wrong results.	The sample is too high concentrated and exceeds the specified absorbance range. Maximum absorbance is 330 A (10 mm path) for nanovolume and 2.65 A for cuvette measurements. Dilute sample and measure again.
Maximum level exceeded.	This message is shown when with an Add + button too many e.g. dyes, wavelength etc. are added. Up to 20 e.g. dyes, wavelength etc. are possible to add.
Measurement currently in progress. Please close method on device.	On the NanoPhotometer is an open measurement session and the message is shown on a control device. Please close method on NanoPhotometer.
No printer present.	Printer connection is lost. Please check printer connection. If it is still not working go back to home screen reconnect the printer and wait for 30 seconds. If the message still appears please create a log file and send it to the Implen customer support.
Sample concentration too low for 0.3µl - change volume and parameter settings to 1µl	The parameter setting volume 0.3 µl measures only the 0.07 mm path length (dilution 140). The sample is too low concentrated use 1 µl and change the volume setting to 1 - 2 µl. Minimum concentrations are for the 0.07 mm path length for dsDNA 420 ng/µl and for BSA 12.6 mg/ml.

### IMPORTANT ALERT MESSAGES:

Cuvette port blocked - please check.	The automatic sliding cover of the NanoPhotometer NP80 could not close properly. Please check if the cuvette port is empty. If message still appears please contact the Implen customer support.
ERROR: Lid open - please close lid.	Lid is open during initialization test. Close lid and confirm with OK
Firmware not found - NPOS.bin must be in root folder of USB flash drive	Make sure that the sample name is NPOS.bin without any additions or extensions like from multiple downloads, otherwise rename file to be NPOS.bin exactly. NPOS.bin file must be in the root folder of the USB flash drive.
Initialization failed – problem with optical path. Please contact Implen customer support.	Check if the cuvette port is empty and/or clean the nanovolume pedestal and the mirror in the lead with 70% EtOH and distilled water and reboot the NanoPhotometer. If the message still appears please create a log file and send it to the Implen customer support.

Initialization failed 51/52/53/54/55. Please contact Implen customer support.	There might be some hardware failure. Please create a log file and send it to the Implen customer support.
Initialization failed. Please contact Implen customer support.	Clean the nanovolume pedestal and the mirror in the lead with 70% EtOH and distilled water and reboot the NanoPhotometer. If the message still appears please create a log file and send it to the Implen customer support.
Light path blocked – please remove cuvettes from port.	Make sure that the cuvette port is empty during initialization test. Restart the NanoPhotometer with empty cuvette port. If the message still appears please create a log file and send it to the Implen customer support.
Measurement head not clean - please wipe pedestal and lid.	Clean the nanovolume pedestal and the mirror in the lead with 70% EtOH and distilled water and reboot the NanoPhotometer. If the message still appears please create a log file and send it to the Implen customer support.
No connection possible	Connection is lost between control devices and NanoPhotometer. Check the connection. If the message still appears please create a log file and send it to the Implen customer support.
Server Not Available	Internal software communication problem during an open measurement session. Reboot the NanoPhotometer. If the message still appears please create a log file and send it to the Implen customer support.
Server Side Failure	Internal software communication problem during an open measurement session. Reboot the NanoPhotometer. If the message still appears please create a log file and send it to the Implen customer support.
Software update failed. Please contact Implen customer support.	Make sure that the sample name is NPOS.bin without any additions or extensions like from multiple downloads, otherwise rename file to be NPOS.bin exactly. NPOS.bin file must be in the root folder of the USB flash drive. Try the update a second time, if it still fails please create a log file and send it to the Implen customer support.
Stored method/result was created for unavailable mode NanoVolume/Cuvette/Submicroliter Cell!	Message appears if a stored results or stored method made on another NanoPhotometer type is opened with an unavailable mode like Cuvette or NanoVolume e.g. data or methods made in cuvette mode and opened on a N60 (NanoVolume only).
Cuvette port blocked - please check.	The automatic sliding cover of the NanoPhotometer NP80 could not close properly. Please check if the cuvette port is empty. If message still appears please contact the Implen customer support.

## 9. WARRANTY

Implen guarantees that the product supplied has been thoroughly tested to ensure that it meets its published specification. The warranty included in the conditions of supply is valid for 12 months and is valid only if the product has been used according to the instructions supplied. Implen or your supplier can accept no liability of loss or damage arising from the faulty or incorrect use of this product

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