ENGLISH: IMPLEN OD600® INSTALLATION

CZECH: INSTALACE IMPLEN OD600®

DUTCH: IMPLEN OD600® INSTALLATIE

FRENCH: IMPLEN OD600® INSTALLATION

GERMAN: IMPLEN OD600® INSTALLATION

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IMPLEN OD600® INSTALLATION

Safety Information

Before commencing installation, please take time to familiarize yourself with warning labels and symbols on your instrument and their meaning. These are to inform you where potential danger exists or particular caution is required.

=== Direct current

Do not open the instrument as this can expose the operator to electrical power and delicate optical setup or damage the instrument.

Do not use damaged power cords, accessories, and other peripherals with your Implen OD600®. Use only the delivered and specified power supply/charger.

Do not expose the Implen OD600® to strong magnetic, electrical fields, water or any type of liquid as heavy rain or moisture.

Do not put the instrument into fire, as it may swell or explode (battery). Do not store at or use near any type of heat source, especially temperatures above 50°C.

Do not leave your Implen OD600® on your lap or near any part of your body to prevent discomfort or injury from heat exposure.

Do not place objects on top of the Implen OD600®.

The Implen OD600® has to be switched off during transportation.

 The symbol on the product, or on the documents accompanying the product, indicates that this appliance may not be treated as household waste. Instead it shall be handed over to the applicable collection point for the recycling of electrical and electronic equipment. Disposal must be carried out in accordance with local environmental regulations for waste disposal.

Unpacking and Positioning

Check the contents of the package against the delivery note. If any shortages are discovered, inform your supplier immediately.

Inspect the instrument for any signs of damage caused in transit. If any damage is discovered, inform your supplier immediately.

Ensure your proposed installation site conforms to the environmental conditions for safe operation: indoor use or dry environment.

The Implen OD600® is designed and developed for use in a controlled, secure environment. This is described as a laboratory environment with restricted physical access within a company. The Implen OD600® may only be accessed by qualified users and may only be operated within a local company network. If a secure environment is not ensured, the Implen OD600® does not meet the following requirements of the Radio Equipment Directive (RED DA): ACM, AUM, SUM, SSM, SCM, CCK, GEC.

Note: Do not expose your Implen OD600® to liquids, rain, moisture or dusty environments.

Operational temperature range: 15 – 40°C

Storage temperature range 5 – 50°C

Transport temperature range: -20°C to 50°C

If the instrument is subjected to extreme temperature changes, it may be necessary to allow the instrument to equilibrate and to turn the instrument off and then on again once thermal equilibrium has been established (~2 – 3 hours).

Maximum relative humidity (non-condensing) of 80% and up to 31°C decreasing linearly to 50% at 40°C.

The instrument must be placed on a stable, level surface that can support 0.6 kg and so that air can circulate freely around the instrument. Ensure while powered on that no materials reduce air circulation.

The equipment should be positioned such that in the event of an emergency the main plug can be easily located and removed.

Carry the instrument always by holding the main corpus of the instrument.

The equipment must be connected to power with the USB-C charger – output: 20W max., input: AC 100-240V, 50/60Hz, 0.7A. The power supply/cord supplied by Implen is a USB-C adapter with 5V. Please charge the battery for at least for 3 hours prior to the first use.

Turn the instrument on using the power button on the touch-panel after it has been plugged in. The instrument will perform a series of self-diagnostic checks.

Please read the complete user manual prior to use.

Please contact original supplier immediately if technical or sample handling difficulties are experienced.

Note: If this equipment is used in a manner not specified or in environmental condition not suitable for safe operation, the protection provided by the equipment may be impaired and the instrument warranty voided.

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